



400 MHz US /R

● User Manual



This Manual was written by
Peter van Bommel

Copyright by Bruker BioSpin AG

All rights reserved. No part of this publication may be reproduced, stored in a retrieval system, or transmitted, in any form, or by any means without the prior consent of the publisher. Product names used are trademarks or registered trademarks of their respective holders.

© November 2010, Bruker BioSpin AG
Fällanden, Switzerland
ZTKS0113 / Z31821 / Rev.: 02

0	Contact	9
1	Introduction.....	11
2	Safety	15
3	Transportation	31
4	Assembling	37
5	Operation.....	39
6	Troubleshooting	41
7	Maintenance.....	57
8	Disassembling	61
A	Appendix	63



0	Contact	9
1	Introduction.....	11
1.1	General Information.....	11
1.1.1	Limitation of Liability	11
1.1.2	Customer Service.....	11
1.1.3	Warranty.....	11
1.1.4	Copyright.....	11
1.2	General View.....	12
2	Safety.....	15
2.1	Persons in charge	15
2.2	Customer Responsibilities.....	16
2.3	Key Words.....	17
2.4	Residual Risks.....	18
2.4.1	Persons	18
2.4.2	Intended Use	18
2.4.3	Safety Devices	19
2.4.4	Spare Parts	19
2.4.5	Signs and Labels.....	20
2.4.6	Technical Risks	20
2.4.7	Cautions	25
2.5	Personal Protective Equipment.....	26
2.6	Description of Signs and Labels.....	27
2.7	Safety Devices	28
2.8	Behavior in Danger and Emergency Situations.....	29
2.9	Fire Department Notification.....	29
3	Transportation	31
3.1	Safety	31
3.2	Packaging.....	31
3.2.1	Disposal.....	32
3.3	Transport Inspection.....	32
3.4	Transportation by Fork Lift / Pallet Jack	33
3.5	Transportation with a Crane	34
3.6	Storing	35

4	Assembling	37
4.1	Safety	37
5	Operation.....	39
5.1	Safety	39
5.2	Set into Operation.....	40
6	Troubleshooting	41
6.1	Safety	41
6.2	Problem	42
6.2.1	During Transportation.....	42
6.2.2	During Operation of the Magnet Stand.....	43
6.2.3	During Standard Operation.....	45
6.2.4	During Operation of the Cryogenic Refrigerator	48
6.3	Troubleshooting Work	53
6.3.1	After a quench	53
6.3.2	Procedure in case of an alarm signal	54
6.3.3	Procedure in case of Cryogenic Refrigerator failure.....	55
7	Maintenance.....	57
7.1	Safety	57
7.2	Cleaning	58
7.3	Maintenance Timetable	58
7.4	Maintenance Work at the Cryogenic Refrigerator	60
8	Disassembling	61
8.1	Safety	61
A	Appendix	63
A.1	Warning Signs	65
A.2	Figures.....	67
A.3	Tables.....	69
A.4	Glossary / Abbreviations.....	71
A.5	Technical Data MS 400'54 USR.....	73
A.6	Index.....	101





0 Contact

Manufacturer:

Bruker BioSpin AG

Industriestrasse 26

CH-8117 Faellanden

Sitserland

Phone: +41-44-825-91-11

Fax: +41-44-825-96-96

<http://www.bruker.com>

Please refer to the Model No., Serial No. and Internal Order No. in all correspondence regarding the NMR system or components thereof.

1 Introduction

1.1 General Information

This manual contains important information about the handling of the supplied magnet system as a part of the magnet system used for NMR analyses. The compliance with all safety and handling instructions, the applicable local accident prevention and general safety regulations are necessary for safe work.

The manual is part of the product. It must be kept in the immediate vicinity of the magnet system and unimpeded access must be ensured at any time.

Read the manual carefully before handling the magnet system or its components.

1.1.1 Limitation of Liability

The information in this manual will take into account the current state of the technology.

The manufacturer assumes no liability for damages resulting from:

- non-compliance with the instructions and all applicable documentation,
- use for purposes not intended,
- not sufficiently approved persons,
- arbitrary changes or modifications and
- use of unauthorized spare parts or accessories.

1.1.2 Customer Service

Technical support is provided by Bruker service via telephone or e-mail. For contact information see page 9 of this document.


1.1.3 Warranty

The warranty terms can be found in the sales documents of the magnet system and in the Terms and Conditions of Bruker BioSpin AG.

1.1.4 Copyright

No part of this publication may be reproduced, stored in a retrieval system, or transmitted, in any form or by any means without the prior consent of the publisher. Product names used are trademarks or registered trademarks of their respective holders.

1.2 General View

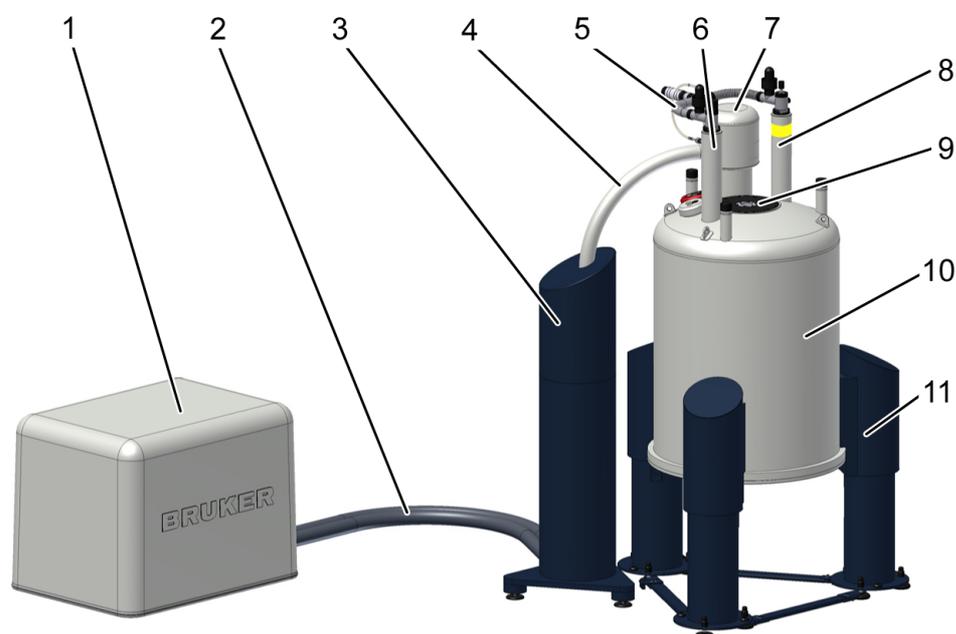

- 1 Cryogenic Refrigerator Compressor (noise protection cover optionally)
- 2 Cryogenic Refrigerator Flex Lines
- 3 Rotary Valve (RV) covered by the Rotary Valve Column (RVC)
- 4 Connection Line
- 5 Helium flow system
- 6 Current lead turret
- 7 Cryogenic Refrigerator Cold Head
- 8 Helium fill-in turret with helium fill-in port

- 9 RT bore
- 10 RT vessel
- 11 Magnet Stand

**HELIUM
FILL-IN PORT**

CAUTION: Use only this port for helium refill!
Don't leave the helium ports open
for more than 5 seconds!
(cf. magnet manual for further information)

Figure 1.1: General view Cryostat of the Magnet System

The heart of the NMR magnet system is a superconducting magnet located inside the helium vessel, which is filled with liquid helium. The helium vessel is surrounded by a radiation shield (RS) cooled by a Cryogenic Refrigerator. The outer casing, the room temperature (RT) vessel (10) contains the helium vessel and the radiation shield. The vacuum inside the RT vessel reduces thermal conduction.

The RT bore (9) allows the access to the magnetic centre of the NMR magnet system. RT vessel, helium vessel, radiation shield, helium turrets, helium flow systems and the RT bore together build the cryostat of the NMR magnet system.

The helium turrets (5, 8) build the interface of the helium vessel to the outside. The current lead turret (5) builds the interface for energizing the magnet coil and for diagnostic. The helium fill-in turret (8) with access to the siphon and connecting plug for the helium level sensor is used to fill the helium vessel. It is marked with a yellow label. The helium flow system (6) connects the two helium turrets.

The Cryogenic Refrigerator is a closed loop helium expansion cycle. It consists of a compressor (1), two flex lines (2), a remote rotary valve inside the rotary valve column (RVC; 3), a cold head (7) and a connection line (4) between rotary valve and cold head. The rotary valve, flex lines, connection line and cold head are covered with an applicable noise protection cover. For the compressor a noise protection is available optionally. Depending on the customers site restrictions several opportunities for the additional equipment of the Cryogenic Refrigerator are possible. Refer to the order subscription and to the supplied manuals of the supplied equipment.

The NMR magnet system is mounted on a magnet stand (8). The isolators in the magnet stand absorb floor vibrations. Different heights and isolators are available optionally.

2 Safety

The supplied cryostat of the magnet system was designed and manufactured according best available technical knowledge and practice, archived in over 50 years of experience of Bruker BioSpin AG. International standards for quality and approval recommended for cryostats of superconducting magnets were certified.

Nevertheless non-compliance with the following instructions and safety advices may cause serious hazards and property damage.

2.1 Persons in charge

In this manual the following qualifications are referenced:

Bruker Service

As a result of professional training, experience and knowledge of applicable regulations they are approved to perform work on magnet systems. They are approved to self-identify dangerous situations and to take applicable measures.

Approved Persons of the Customer

As a result of professional training, experience and knowledge of applicable regulations and instruction by the customer, this group of people is approved to perform the duties assigned to them. They are approved to recognize dangerous situations and to take applicable measures.

2.2 Customer Responsibilities

The customer of the magnet system must obey the security advice and the rules for safety, accident prevention and environmental protection correctly for the magnet systems. Furthermore, the customer is responsible for keeping the magnet system in correct technical condition.

In particular:

- The customer must determine additional dangers resulting from the working conditions at the site of the magnet system and provide applicable safety measures.
- The customer must ensure that the site plan meets the specified conditions for operating the magnet system and supplied site planning document.
- The customer has to mark the danger area around the magnet system and post the corresponding instruction plates.
- The customer has to inform the local fire brigade about the special risks of the magnet system and how to react in the event of an incident.
- The customer must clearly define the responsibilities for operation and maintenance.
- The customer must ensure that all employees working with the magnet system have read and understood the manual.
- The customer has to provide the necessary personal protective equipment for his employees.
- The customer has to instruct his employees at regular intervals on hazards and safety measures.
- The customer has to instruct other persons not working on the magnet system but carrying out work in the same room, for instance cleaning staff.
- The customer has to consider the specific items of this N₂-free cryostat equipped with a Cryogenic Refrigerator. The customer is responsible for obeying the advice given in this manual. In case the Cryogenic Refrigerator is not running correctly his immediate reaction is mandatory. In case of an unexpected alarm his immediate response is mandatory. For further instruction refer to chapter "**Troubleshooting**" on page 6-41.
- The customer has to obey the specified maintenance intervals given in this manual (see chapter "**Maintenance Timetable**" on page 7-58).

2.3 Key Words



DANGER

Indicates a hazardous situation which, if not prevented, will result in death or serious injury.



WARNING

Indicates a hazardous situation which, if not prevented, could result in death or serious injury.



CAUTION

Indicates a hazardous situation which, if not prevented, may result in minor or moderate injury.

NOTICE

Hazard, which could result in property damage.



Information and links for efficient and trouble-free handling and operation.

2.4 Residual Risks

In the following section, the residual risks from the risk analysis according ISO 14971 are summarized. To prevent health hazards and hazardous situations obey all safety instructions and warnings in the manual.

2.4.1 Persons

WARNING



Risk of injury and property damage from not approved persons.

Incorrect handling of the magnet system from not approved persons may result in significant bodily injury and property damage.

Thus:

- Work must only be carried out by approved persons with applicable qualifications. The necessary qualifications are specified in the beginning of the relevant chapter.
- In case of doubt, contact Bruker Service. Contact information see page 9 of this document.

2.4.2 Intended Use

WARNING



Risk of damage to life and limb by incorrect use of the magnet system.

Incorrect use of the magnet system can lead to life-threatening situations and destruction of the magnet system.

Thus:

- Only use the magnet system as intended.
- Do not change the magnet system.
- Do not use the magnet system to demonstrate strong magnetic fields.
- Do not exceed specified values for operating the magnet system.

Damage claims from damages caused by other than the intended use of the magnet system are excluded and the customer is held liable.

2.4.3 Safety Devices

WARNING



Risk of damage to life and limb due to not sufficient safety devices.

Several safety devices ensure safe operation of the magnet system. They must always be in correct working condition.

Thus:

- Do not block safety devices.
- Do not remove safety devices.
- Check the operational reliability of the safety devices before working on the magnet system.

2.4.4 Spare Parts

WARNING



Risk of injury and property damage from using incorrect or defective spare parts and accessories.

Incorrect or defective spare parts can cause serious injuries. They may cause damaging, malfunctioning and the destruction of the magnet system.

Thus:

- Use only original equipment manufacturer spare parts.
- Use only original equipment manufacturer accessories.

2.4.5 Signs and Labels



WARNING

Risk of damage to persons and property due to not readable signs and labels.

Signs and labels with advice may become not readable.

Thus:

- Maintain signs and labels in a readable state.
- Replace damaged or not readable signs and labels immediately. New signs and labels can be obtained from Bruker Service.

2.4.6 Technical Risks

Magnetic Field



WARNING

Risk of damage to life and limb due to high magnetic fields.

A magnetic field of more than 0.5 mT (5 Gauss) is life-threatening for people with pacemakers or metal implants. Ferromagnetic tools in the magnetic field are significantly hazardous. Disks and electronic devices may be damaged. Duration of exposure (8 h/day) above the limit of 200 mT can cause damage to health.

Thus:

- Mark the magnetic field of more than 0.5 mT (5 Gauss) before start up.
- The workplace must be outside the 0.5 mT area.
- Keep people with pacemakers and metal implants away from the identified area.
- Keep disks, credit cards and electronic devices away from the identified area.
- Do not use ferromagnetic tools or items within the identified area.
- Only use non-ferromagnetic transportation dewars for the cryogenic agents.
- Only use non-ferromagnetic ladders or steps.

Cryogenic Agents

WARNING



Risk of damage to life and limb due to cryogenic agents.

Risk of damage to life and limb due to not correct handling of liquid cryogenic agents and their dewars. Within the transition from liquid to gas, helium and nitrogen expands their volume, causing closed vessels to burst.

The evaporating cryogenic agents will displace the air. In case of not sufficient ventilation this may result in death by suffocation. Helium displaces oxygen in the upper part of the room, nitrogen displaces oxygen in the lower parts of the room.

Liquid and gaseous cryogenic agents are extremely cold. Contact with liquid or gaseous cryogenic agents will lead to cold burns.

Thus:

- Only use cryogenic agents in well ventilated rooms. In case of doubt ask Bruker Service.
- Wear an oxygen monitor on the body during service and maintenance work.
- Prevent any skin contact with liquid or gaseous cryogenic agents.

Electricity

WARNING



Risk of damage to life and limb due to electricity.

Risk of damage to life and limb due to contact with electrical lines and damaged insulation!

Thus:

- Work on electrical equipment must be done by an approved electrical technician.
- Keep moisture away from electrical lines to prevent short-circuits.
- Check the magnet system accidental ground before start up.
- Switch power off before working on the power supply or further equipment.

Quench



WARNING

Risk of suffocation during a quench of the magnet system.

A quench is the very fast de-energizing of the magnet by loss of its superconductivity. The stored magnetic energy is converted into heat and thus large quantities of helium evaporate. The evaporating helium will displace oxygen. In case of not sufficient ventilation this may result in death by suffocation.

Thus

- The magnet system site must be well ventilated. In case of doubt contact Bruker Service.
- The evaporating gas may resemble smoke. Never pour water on the magnet system.

Gas under Pressure



WARNING

Risk of injury due to gas under pressure inside the cryostat and further equipment.

The helium vessel of the cryostat may get sealed off due to ice formation inside the helium turrets in case of non-compliance with the instruction given in this manual. This may lead to overpressure and damage of the helium vessel.

Manipulations of components with gas under pressure may lead to injury and property damage.

During cold head warm up the helium in the Cryogenic Refrigerator circuit gas expands. If the cryo cooler circuit is open, the pressure in the system increases. The overpressure will be released through the security valves on cold head and rotary valve. The expelled gas can cause severe cold burns.

Thus:

- In case of icing inside the helium turrets contact Bruker Service immediately.
- Release the pressure before working on components with gases under pressure.
- Do not seal cryogenic agent vessels of the magnet system or the transport dewars.
- Do not connect high pressure transport dewars for liquid cryogenic agents to the magnet system. Completely eliminate the high pressure from the transport dewars before connecting and transferring.
- Keep the Cryogenic Refrigerator circuit closed at any time. Overpressure can release in the buffers of the compressor, the rotary valve and the cold head.

Low Temperatures

WARNING



Risk of injury due to low temperatures of liquids and metal parts.

Physical contact with extremely cold liquids and metal parts may cause serious injuries. Contact with the skin may cause cold burns. Contact with the eyes may cause blindness.

Thus:

- Always wear protective goggles, protective gloves and protective clothes while handling with liquid cryogenic agents or metal parts in contact with cryogenic agents.
- Protect temperature sensitive components such as O-rings from contact with liquid cryogenic agents.

Spontaneous Ignition and Explosion

WARNING



Risk of injury from spontaneous ignition and explosion caused by liquid oxygen.

Pure oxygen condenses on extremely cold metal pieces. Together with oil it may ignite spontaneously. In case of fire the pure oxygen may cause an explosion.

Thus:

- Do not smoke near the magnet system.
- Do not use open flames near the magnet system.
- Keep the environment around the magnet system clean.
- Do not leave oily rags near the magnet system.

Risk of Slippage

WARNING



Risk of injury from slippage.

The accumulation of condensed water on the floor and ladders causes slippery surfaces.

Thus:

- Always wear safety shoes with an anti slip sole.
- Be careful using ladders.
- Clean floor and ladders regularly.

Risk of Tilting



WARNING

Risk of injury due tilting of the magnet system.

The magnet system is very sensitive to lateral forces. It may tilt.

Thus:

- Do not climb onto the magnet system.
- Do not lean items against the magnet system.
- Do not lean against the magnet system.
- Do not move the magnet system on your own.

Risk of Heavy Weights



WARNING

Risk of damage to life and limb caused from heavy moving weights.

Lifting heavy weights is life-threatening due to falling or moving parts.

Thus:

- Do not stay or work under a lifted magnet system.
- All lifting equipment used must be applicable to carry the weight (see Appendix, Technical Data).
- Do not use damaged lifting equipment.
- Do not use lifting equipment without updated check tag.
- Lifting only with approved qualification.
- Obey ergonomic guidelines while lifting heavy parts.
- Protect parts against falling.

Risk of Hot Surfaces



WARNING

Risk of injury from contact with hot surfaces.

Surfaces of the Cryogenic Refrigerator parts may be hot. Skin contact with these surfaces can cause serious injuries.

Thus:

- Any work at the Cryogenic Refrigerator parts must only be carried out by Bruker Service or approved persons.
- Always wear protective gloves.

2.4.7 Cautions

CAUTION

Risk of injury and property damage due to incorrect transportation.

The box may tilt, movement may get out of control. Thus persons may get injured and the cryostat or further equipment may be damaged.

Thus:

- Do not move arbitrarily.
- Be careful while unloading and moving the boxes.
- Pay attention to any symbols on the boxes.
- Move the box in an upright position.
- Do not tilt the box.
- Prevent crossing thresholds, even if they are only a few millimeters high.
- Clean the transportation way before transporting the box.
- Unpack shortly before assembling.
- Only use the attachment points provided.
- Transportation only with transportation locks attached.
- The cryostat or further equipment must be protected from rain and other bad weather conditions during transportation.
- Pay attention to sharp edges and spikes of boxes and parts by using protective gloves while moving.



2.5 Personal Protective Equipment

The personal protective equipment must be worn at any time while working on the magnet system and to avoid health hazards.



Protective Goggles

Used to protect the eyes from injury due to flying cold liquids and parts.



Protective Gloves

Used to protect the hands from injury caused by contact with extremely cold liquids or surfaces and for protection from injury caused by rough edges.



Protective Clothes

Used to protect the body from injury caused by contact with extremely cold liquids or surfaces and for protection from wounds.



Safety Shoes

Used to protect the feet from injury from falling of heavy objects. An anti-slip sole protect from injury caused by slipping and falling on slippery floor and steps. Only use safety shoes with non-ferromagnetic toe caps.



Portable Oxygen Monitor and Alarm

Used to warn against low oxygen concentrations in surrounding air.

2.6 Description of Signs and Labels

Signs and labels are always related to their immediately vicinity. The following signs and labels are found on the magnet system and in the vicinity.



Prohibition sign: No person with pacemakers!

People with pacemakers are endangered in the identified area of 0.5 mT (5 Gauss) and are not allowed to enter these areas.



Prohibition sign: No person with implants!

People with metallic implants are endangered in the identified area of 0.5 mT (5 Gauss) and are not allowed to enter these areas.



Prohibition sign: No watches or electronic devices!

Watches and electronic devices may be damaged in the identified area of 0.5 mT (5 Gauss).



Prohibition sign: No credit cards or other magnetic memory!

Credit cards and magnetic memory may be damaged in the identified area of 0.5 mT (5 Gauss).



Prohibition sign: Do not touch! Do not block!

Do not touch or block identified area.



Hazard warning sign: Strong magnetic field!

- No magnetic memory.
- No jewelry.
- No metallic items.



Emergency exit!

- Always keep the emergency exit clear.
- Follow the arrows if necessary.
- Doors must be pushed open in escape direction.

2.7 Safety Devices

The cryostat of the supplied magnet system is equipped with the following safety devices:

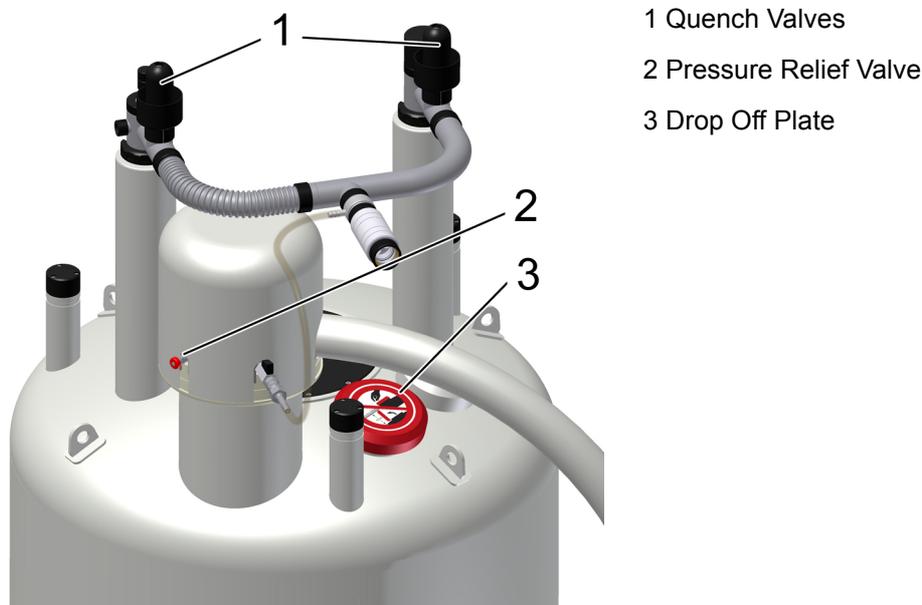


Figure 2.1: Safety Devices of the Cryostat

Quench Valve

The quench valves (1) are the safety devices of the helium vessel. They open with a defined pressure. In case of an accidental overpressure in the helium vessel the quench valves will release the pressure smoothly.

Pressure Relief Valve

The pressure relief valve (2) is the safety device of the cold head turret. It opens with a defined pressure. In case of an accidental overpressure in the cold head turret the pressure relief valve will release the pressure smoothly.

Drop Off Plate

The drop off plate (3) is a safety device of the RT vessel. If the vacuum breaks, the drop off plate will open. In case of an accidental overpressure in the RT vessel the drop off plate will release the pressure smoothly.

Cryogenic Refrigerator Parts

For information about the safety devices of the Cryogenic Refrigerator parts refer to the supplied separate manual.

2.8 Behavior in Danger and Emergency Situations

Preparations

- Keep the emergency exits free at all times.
- Prepare and maintain an up-to-date list of emergency telephone numbers in the magnet system area.

In Case of Emergency

- Leave the danger zone immediately.
- Check for sufficient ventilation in the room before entering, especially if people are showing symptoms of suffocation.
- Rescue persons from the danger zone.
- Provide medical attention for people with symptoms of suffocation.
- Start first aid immediately.
- Call the responsible contact.
- Call for medical assistance.
- Call the fire department.

First Aid for Cold Burns

- Help the injured persons to lie down comfortably in a warm room.
- Loosen all clothing which could prevent blood circulation in the injured area.
- Pour large quantities of warm water over the affected parts.
- Cover the wound with dry and sterile gauze.
- In case of contact of liquid cryogenic agents with the eyes rinse thoroughly with clean water.
- Call for medical assistance.

2.9 Fire Department Notification

- Inform the fire department about the potential risks of a magnet system, i.e. danger of metallic rescue equipment close to the magnet system.
- Laboratory windows which are accessible during an emergency should be clearly identified with warning signs, visible from the outside.
- Inform the fire department about the characteristics of a quench to prevent confusion with smoke.
- Never pour water over the magnet system during a quench!

3 Transportation

3.1 Safety

The transportation is carried out by Bruker Service or approved persons. However, it may happen that other persons have to receive the delivery of the shipment boxes. In this case it is essential to obey the instructions in this chapter and to inform these persons before.



WARNING

Heavy Weights



CAUTION

Incorrect Transportation

3.2 Packaging



Figure 3.1: Packaging (without surrounding panels)

The cryostat is supplied in a wooden box on a pallet. It is secured inside with straps against tilting and moving.

The two flex lines of the Cryogenic Refrigerator are wrapped around the cryostat, fixed with the same straps.

Accessories such as the flow systems, level sensors and bore tubes are in the side compartment of the box.

The Cryogenic Refrigerator parts are supplied in a wooden box on a pallet.

The Magnet Stand is supplied in a wooden box on a pallet.

3.2.1 Disposal

Keep the original boxes for future transportation.

If no further transportation is planned, dispose of the boxes according to environmentally friendly regulations.

3.3 Transport Inspection

Investigate the delivery with regard to visible damage and completeness of delivery.

Delivery list

- Cryostat box
- Cryogenic Refrigerator box
- Magnet Stand box

Transport control systems

The shipping and handling monitors ("Shock Watch", "Tilt Watch") on the boxes shows, if the boxes were kicked or tilted during transportation.

Checks

Shock Watch: Follow instructions on the label.

Tilt Watch: Follow instructions on the label.

In case of damage

- Accept delivery with reservation.
- Make a documentation of all observable damage and add it to the transport documents.
- Start complaint process.
- Contact Bruker Service before installation.



The claim for damage expires after the fixed period

Thus:

Report damages to Bruker Service immediately after detection of damage. For contact information see page 9 of this document.

3.4 Transportation by Fork Lift / Pallet Jack

Recommended for transporting the boxes to the installation site.

Persons in charge: Approved forklift / pallet jack operator

Precondition: The fork lift/pallet jack must be approved for the transportation weight (see chapter Appendix, Technical Data).

Transport



1. Check the route of transport for the minimal height and width.
2. Check sufficient floor capacity on the route of transport. In case of doubt ask a stress analyst.
3. Check sufficient carrying capacity while using an elevator.
4. Position the forks between the bars of the box as shown in the figure. Make sure the side towards the operator is the one with the labels on it.
5. Make sure the forks of the fork lift are longer than the box and projects out of the back of the box as shown in the figure below.



6. Now lift the fork and move the box to the site.

Figure 3.2: Transportation by forklift


3.5 Transportation with a Crane

Recommended for lifting the cryostat out of the box.

Persons in charge: Approved crane operator

Precondition: The crane must be approved for the transportation weight (see chapter Appendix, Technical Data).

Attachment Points

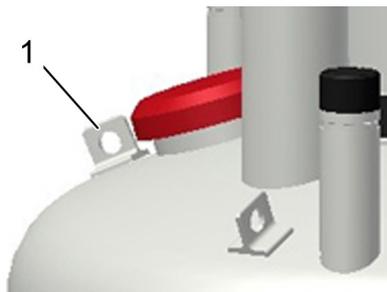

1. Exclusively use the marked eyelets (1) as attachment points for the lifting equipment.
2. Use **all** eyelets for the lifting equipment.
3. Follow the instructions on the label on the top of the cryostat. This label gives important information about correct attachment and transportation.
4. Check for correct fastening of the lifting equipment before lifting the cryostat.
5. Make sure that any movement of the crane is as slowly as possible to avoid any damage due to acceleration.
6. Check for correct leveling of the cryostat while hanging on the crane.

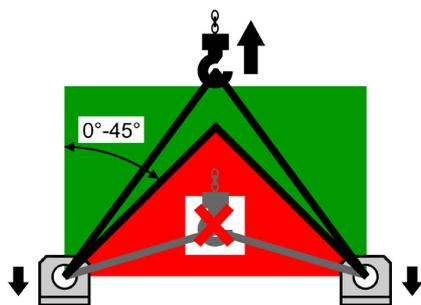


Figure 3.3: Attachment points for lifting equipment

3.6 Storing

If it is necessary to store the cryostat and accessories before installation obey the following instructions:

- Store the boxes in a closed, dry and dust-free room.
- Store the boxes upright.
- Do not tilt the boxes.
- Do not unpack the supplied boxes.
- Prevent mechanical vibrations to the boxes.
- Storage temperature: 5 - 40°C.
- Storage humidity: less than 50%@23°C.

4 Assembling

4.1 Safety

Persons in charge: Bruker Service only

5 Operation

5.1 Safety

WARNING



Gas under Pressure
Cryogenic Agents
Electricity

Persons in charge:

Bruker Service, approved persons of the customer

5.2 Set into Operation

Mount the further equipment of the supplied magnet system respecting their manuals.



Figure 5.1: Start the Magnet Stand

Set the Magnet Stand into operation by switching the pneumatic controller to UP position (1).



Figure 5.2: Stop the Magnet Stand

For any work at the magnet system like maintenance or refill of helium stop the Magnet Stand by switching the pneumatic controller to DOWN position (1).

6 Troubleshooting

Troubleshooting must be performed only with approved qualification.

In case of doubts or problems not specified in the following list contact Bruker Service immediately. For contact information see page 9 of this manual.

6.1 Safety

WARNING



Magnetic Fields

Gas under Pressure

Cryogenic Agents

Electricity

Spontaneous Ignition and Explosion

Persons

Bruker Service, approved persons of the customer

6.2 Problem

6.2.1 During Transportation

Indicator	Possible reason	Solution	By
Tilt Watch / Shock Watch activated.	Careless transportation.	<ol style="list-style-type: none">1. Accept delivery with reservation.2. Remark the extent of damage in the transport documents.3. Start complaint process.	Approved persons of the customer
Visible damage.	Careless transportation.	<ol style="list-style-type: none">1. Accept delivery with reservation.2. Remark the extent of damage in the transport documents.3. Start complaint process.	Approved persons of the customer

6.2.2 During Operation of the Magnet Stand

In case of doubt contact Bruker Service and refer to the supplied manual of the Magnet Stand

Indicator	Possible reason	Solution	By
The NMR spectrum shows massive disturbances.	The pneumatic controller is in DOWN position.	Switch pneumatic controller to UP position.	Approved persons of the customer
	Magnet system has direct mechanical contact with the floor via accessories.	Identify and eliminate contact point. Arrange lines in loose S- or U shapes	Approved persons of the customer
		If the problem is still not solved, contact Bruker Service.	Approved persons of the customer
	Magnet system has physical contact to the magnet stand.	Check leveling; adjust if necessary.	Bruker Service
	Piston of the isolator is not centric or touches its casing.	Align magnet stand.	Bruker Service
	Floor vibrations in horizontal and vertical direction.	Replace isolators with pneumatic isolators with horizontal isolation.	Bruker Service

Continued on next page

Continued from page before

Indicator	Possible reason	Solution	By
The isolator of the magnet stand does not reach the operating position.	Pneumatic controller in DOWN position.	Switch pneumatic controller to UP position.	Approved persons of the customer
	The pressure of the gas supply is too low.	Check the pressure of the gas supply. It must be in the range of 5 to 8 bar (70 to 112 psi).	Approved persons of the customer
		If the problem is still not solved, contact Bruker Service.	Approved persons of the customer
	The magnet system is not leveled correctly.	Stop the pneumatic isolators. Check the leveling of the cryostat.	Bruker Service
	Defective leveling valve.	Exchange leveling valve or isolator.	Bruker Service
	Defective membrane of an isolator.	Exchange leveling valve or isolator.	Bruker Service
Magnet system achieves working position jerkily.	Piston is not centric or touching its casing.	Align magnet stand.	Bruker Service
Audible loss of air	Defective membrane of an isolator.	Exchange leveling valve or isolator.	Bruker Service
	Hose connector is defective or loose.	Insert hoses correctly and tighten screws.	Bruker Service
Velocity of lifting or lowering too high.	Wrong adjustment of the restrictor level valve.	Close restrictor of the level valve completely; then open it a half turn.	Bruker Service

6.2.3 During Standard Operation

Indicator	Possible reason	Solution	By
The helium boil off decrease to zero.	The atmospheric pressure is increasing.	Usual behavior. Watch helium boil off daily.	Approved persons of the customer
	The helium flow system is covered with ice.	Contact Bruker Service immediately! Do not try to remove ice of the helium flow system!	Approved persons of the customer
		If the problem is still not solved, contact Bruker Service.	Approved persons of the customer
	⚠ WARNING: Cryogenic Agents Quench		
	The helium flow system or the suspension tubes are blocked with ice.	Blow in warm helium gas carefully through an applicable tube. Do not insert it more than 600mm from the top of the helium turrets.	Bruker Service
The helium boil off is too high.	The helium level meter is permanently on (service mode) or used often.	Switch off helium level sensor. Reduce helium level measurement (during measuring of the helium level an amount of helium boils off due to the heat input of the helium level sensor; for reporting the helium boil off a weekly measurement is sufficient).	Approved persons of the customer
	The atmospheric pressure is decreasing.	Usual behavior. Watch helium boil off weekly.	Approved persons of the customer
If the problem is still not solved, contact Bruker Service.		Approved persons of the customer	

Continued on next page

Continued from page before

Indicator	Possible reason	Solution	By
Continue of: The helium boil off is too high.	Vacuum reduced.	Rebuild vacuum, see chapter "Rebuilding the Vacuum" on page 4-53	Bruker Service
	The radiation baffles are not inserted in the current lead turret.	Insert the radiation baffles into the current lead turret.	Bruker Service
RS temperature shield to high.	Cryogenic Refrigeration not sufficient.	See chapter "During Operation of the Cryogenic Refrigerator" on page 6-48.	Approved persons of the customer
	Alarm default settings not correct.	Check set values (see separate manual of the External Temperature Control Unit).	Approved persons of the customer
		If the problem is still not solved, contact Bruker Service.	Approved persons of the customer
	Vacuum reduced.	Rebuild vacuum, see chapter "Rebuilding the Vacuum" on page 4-53.	Bruker Service
Quench	Loss of superconductivity.	See chapter "After a quench" on page 6-53 Contact Bruker Service immediately!	Approved persons of the customer
Cold spots inside the RT bore.	Alignment not correct.	Contact Bruker Service.	Approved persons of the customer
RT vessel is wet and cold.	Vacuum reduced.	Contact Bruker Service immediately!	Approved persons of the customer
Not correct helium level warning out of supplied NMR software.	Helium level sensor defective.	Contact Bruker Service immediately!	Approved persons of the customer
Helium level at constant level, no change during days.	Helium level sensor defective.	Contact Bruker Service immediately!	Approved persons of the customer

Continued on next page

Continued from page before

Indicator	Possible reason	Solution	By
		⚠ WARNING: Low Temperature	
Not correct helium level warning out of supplied NMR software.	Helium level sensor defective.	Replace helium level sensor (see chapter "Replacement of the Helium Level Sensor" on page 6-110)	Bruker Service
Helium level at constant level, no change during days.	Helium level sensor defective.	Replace helium level sensor (see chapter "Replacement of the Helium Level Sensor" on page 6-110)	Bruker Service

6.2.4 During Operation of the Cryogenic Refrigerator

i In the case of a complete failure of the Cryogenic Refrigerator, the helium contained in the helium vessel will be sufficient to cool the magnet for at least 3 days. If solving the issue takes longer, it is necessary to refill helium at all time (refer to supplied Refill Manual).

i In case of any issue or failure specified in the following table where the intervention of the operator is necessary, make sure the start button of the compressor is switched off and the power supply of the compressor is disconnected.
 After resolving the issue, connect the power supply and push the start button of the compressor.
 If the compressor does not start immediately, press the temperature and the pressure switch for reset and after this the start button on the compressor.
 For problems not specified in this chapter refer to the supplied manual of the Cryogenic Refrigerator.

Indicator	Possible reason	Solution	By
Compressor is not operating.	No power supply or power supply interrupted.	Check the power supply to the compressor and verify that it meets the system requirements.	Approved persons of the customer
	Circuit breaker OFF.	Check that the circuit breaker on the front panel of the compressor is ON. Check for possible causes why the circuit breaker switched off.	Approved persons of the customer
		Check for possible causes why the circuit breaker switched off.	Approved persons of the customer
	Pressure switch tripped	Reset the pressure switch located at the bottom of the front panel of the compressor.	Approved persons of the customer
	Temperature switch tripped	Reset the temperature switch located at the bottom of the front panel of the compressor.	Approved persons of the customer

Continued on next page

Continued from page before

Indicator	Possible reason	Solution	By
Compressor is operating, but no pressure fluctuation visible (needles at the pressure gauges not oscillating).	Motor cord disconnected at the compressor.	Connect the motor cord to the compressor.	Approved persons of the customer
	Motor cord disconnected.	Check the motor cord connection at the coupling in the middle of the motor cord at 10m (skip, if motor cord is one piece).	Approved persons of the customer
		⚠ WARNING: Hot Surfaces	
	Motor cord disconnected at the rotary valve.	<ol style="list-style-type: none"> 1. Dismount the noise protection cover of the rotary valve column. 2. Check the motor cord connection at the rotary valve. 3. Connect motor cord, if necessary. 4. Remount the noise protection cover. 	Approved persons of the customer
		If the problem is still not solved, contact Bruker Service.	Approved persons of the customer
	Aeroquip® connectors of the flex lines not correctly tightened at the compressor.	Tighten the Aeroquip® connectors correctly.	Bruker Service
	High and low pressure helium flex line reversed at the compressor.	Check if the flex lines are mounted correctly at the compressor respecting high and low pressure port.	Bruker Service
Motor cord defective.	Check continuity of all 4 (four) conductors in the motor cord. If not correct, change the motor cord.	Bruker Service	

Continued on next page

Troubleshooting

Continued from page before

Indicator	Possible reason	Solution	By
Continue of: Compressor is operating, but no pressure fluctuation visible (needles at the pressure gauges not oscillating).	Aeroquip® connectors of the flex lines tightened not correctly at the rotary valve.	Tighten the Aeroquip® connectors correctly.	Bruker Service
	High and low pressure helium flex line reversed at the rotary valve.	Check if the flex lines are mounted correctly at the rotary valve respecting high and low pressure port.	Bruker Service
	Rotary valve is defective.	If possible, check if the rotary valve sound has changed since installation. If the typical sound is missing, the rotary valve might be defective and has to be exchanged. Replace the rotary valve.	Bruker Service
Compressor is operating, pressure fluctuation visible at the pressure gauges, extinguish short after start.	Water chiller not running.	Check possible reasons why the water cooler is not running. Start the water cooler.	Approved persons of the customer
	Cooling water flow too low.	Check cooling water supply is sufficient (refer to supplied manual of the Cryogenic Refrigerator).	Approved persons of the customer
	Cooling water flow too low due to icing, blockage, fouling or leak in the cooling water lines.	Search possible reason for the disturbance. Check the cooling water supply is as specified in the supplied manual of the Cryogenic Refrigerator and the manual of the water chiller.	Approved persons of the customer
	Not correct environmental temperature (too high or too low).	Check the environmental temperature is as specified.	Approved persons of the customer
	Not correct water temperature (too high or too low).	Check the cooling water temperature is as specified.	Approved persons of the customer

Continued on next page

Continued from page before

Indicator	Possible reason	Solution	By
Compressor is operating, cryogenic refrigeration not sufficient.		Exclude all previous reasons of this list.	Approved persons of the customer
	Not correct maintenance.	Check periodic maintenance work was made according to schedule (see chapter "Maintenance Timetable" on page 7-58).	Approved persons of the customer
		If the problem is still not solved, contact Bruker Service.	Approved persons of the customer
		⚠ WARNING: Hot Surfaces	
	Connection line between rotary valve and cold head not mounted correctly at the cold head.	Tighten the Aeroquip® connectors correctly. Check if the O-rings of the fitting are in correct position (see chapter "Mounting the Rotary Valve and Flex Lines" on page 4-61).	Bruker Service
	Low helium pressure inside the compressor helium circuit.	Recharge helium of high purity (He 6.0, 99,9999%). Refer to supplied manual of the Cryogenic Refrigerator.	Bruker Service
	Low helium pressure due to small leak inside the helium circuit.	<ol style="list-style-type: none"> 1. Leak detection with leak detector at connectors, flex lines, inside compressor, cold head, rotary valve, connection line. 2. Eliminate leak, if possible. 3. Recharge helium of high purity (He 6.0, 99,9999%). Refer to supplied manual of the Cryogenic Refrigerator. 	Bruker Service
High helium pressure inside the compressor helium circuit.	Vent helium to the set value. Refer to supplied manual of the Cryogenic Refrigerator.	Bruker Service	

Continued on next page

Continued from page before

Indicator	Possible reason	Solution	By
Continue of: Compressor is operating, cryogenic refrigeration not sufficient.	Cold head contaminated.	Remount the cold head. Pump and flush the cold head. See chapter " Procedure in case of Cryogenic Refrigerator failure " on page 6-55	Bruker Service
	Cold head defective.	Replace the cold head. See chapter " Procedure in case of Cryogenic Refrigerator failure " on page 6-55	Bruker Service
	Cryogenic Refrigerator contaminated.	Replace the complete Cryogenic Refrigerator. See chapter and " Mounting the Cryogenic Refrigerator Parts " on page 4-54.	Bruker Service

6.3 Troubleshooting Work

6.3.1 After a quench



Figure 6.1: Quench picture

A quench is the very fast deenergizing of the magnet by loss of its superconductivity. The stored magnetic energy is converted into heat, which promotes rapid evaporation of large quantities of helium. After an appearance of a quench contact Bruker Service immediately.



WARNING

Quench Cryogenic Agents

Quench while in persistent mode:

1. Wait until the quench valves are closed and no helium evaporate out of the quench valves.
2. Wait until there is no helium vapor visible in the room or the ceiling to make sure there is sufficient oxygen in the room.
3. Switch off the alarm at the External Temperature Control Unit.
4. Check the globes in the quench valves for their correct position.
5. Remove probe and shim system to prevent icing from the shim system.
6. Start the refill with liquid helium as soon as possible after the quench (within one hour after the quench; refer to the supplied refill manual).
7. Contact Bruker Service immediately.



If the quench occurs unattended or helium transfer was not possible within one hour after the quench, it is recommended to warm up the system to 90 K.

6.3.2 Procedure in case of an alarm signal

If the temperature of the radiation shield exceeds the set maximum allowed temperature, an alarm will warn the operator (audible from the External Temperature Control Unit, visible from the External Temperature Control Unit and the console via MICS).

Follow the procedure to fix the issue before contacting Bruker Service.

Note:

- Date and time
- Temperature given on the display of the External Temperature Control Unit
- Turn the alarm off (see supplied manual of the External Temperature Control Unit)
- Which errors or warnings appear?
- What are the readings of the pressure gauges, run time counter etc. of the compressor?
- Which parts seems to be ok/not ok?
- How did the failure occur and what happened before failure?

6.3.3 Procedure in case of Cryogenic Refrigerator failure

In case of a failure of the Cryogenic Refrigerator the temperature of the shield T_{RS} will slowly rise to 200 K and the helium boil off to 250 ml/h. This is usual behavior.

In case of failure shorter than 3 days just restart the compressor of the Cryogenic Refrigerator. See remarks at chapter "During Operation of the Cryogenic Refrigerator" on page 6-48 for further information.

After restarting the Cryogenic Refrigerator, at least two days are necessary to recover the previous state of the magnet system. If the temperature and helium boil off does not decrease in the next two days call Bruker Service.

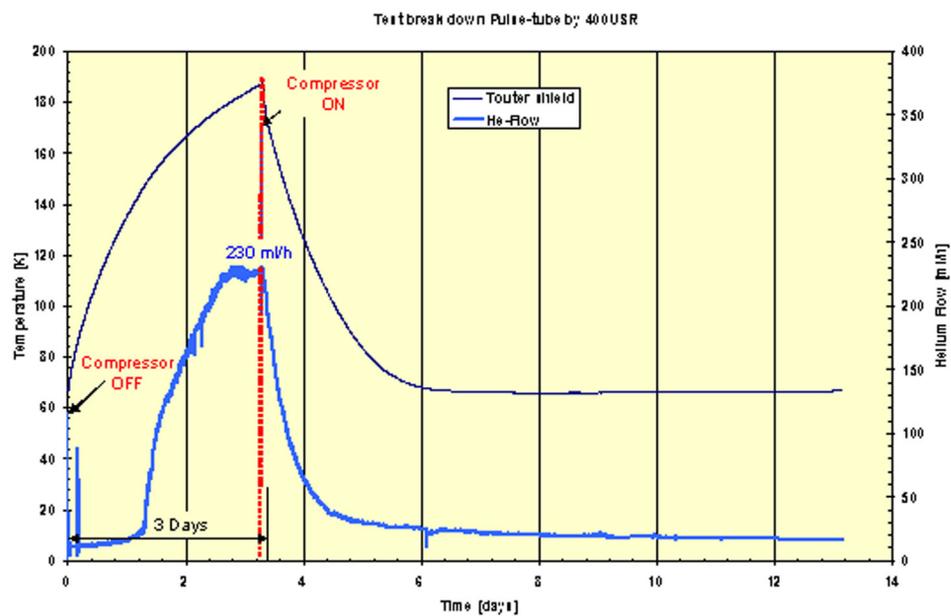


Figure 6.2: Temperature profile during Cryogenic Refrigerator failure

7 Maintenance

Maintenance must be performed only with approved qualification.

In case of doubt contact Bruker Service. For contact information see page 9 of this document.

7.1 Safety

WARNING



Magnetic Fields

Cryogenic Agents

Electricity

Gas under Pressure

Low Temperatures

Spontaneous Ignition and Explosion

Hot Surfaces

Persons in charge

Bruker Service, approved persons of the customer

Personal protective equipment

Protective goggles

Protective gloves

Safety shoes

7.2 Cleaning

Procedure

- Clean the RT vessel of the magnet system, the magnet stand with dry or slightly damp cloth.
- Only use neutral water and detergents.
- Do not use volatile cleaning solvents.

7.3 Maintenance Timetable

- i** If service is required, contact Bruker Service immediately to reduce possible down time due to lack of spare parts or liquid helium supplies.

Interval	Device	Work	By
weekly	Cryostat	Check the helium flow.	Approved persons of the customer
4.000 h (0.5 year)	Cryogenic Refrigerator	<ul style="list-style-type: none"> • Check the values at the high and low pressure gauges on the front panel of the compressor if the mean values are different and if an oscillation of the needles is noticeable. Refer to the supplied manual of the cryogenic refrigerator. Contact Bruker Service if values are different. <p>Water cooled option only:</p> <ul style="list-style-type: none"> • Check the proper operation of the water cooling unit. • Check system pressure, water flow and temperatures. Refer to the manual of the water cooling unit for further information. 	Approved persons of the customer
8.760 h (1 year)	Cryostat	<ul style="list-style-type: none"> • Refill helium. Refer to supplied Refill Manual. • Record the filling session. 	Approved persons of the customer

Table 7.1: Maintenance Timetable - part 1

Interval	Device	Work	By
17.520 hours (2 years)	Cryogenic Refrigerator	<ul style="list-style-type: none"> Replace the rotary valve. Refer to chapter "Replacement of the Rotary Valve and the Cold Head" on page 7-118. Replace the adsorber of the compressor. Refer to the supplied manual of the cryogenic refrigerator. Check the values at the high and low pressure gauges on the front panel of the compressor if the mean values are different and if an oscillation of the needles is noticeable. Refer to the supplied manual of the cryogenic refrigerator. <ul style="list-style-type: none"> - Pressure to high: Vent helium gas. Use the Service Kit. - Pressure to low: Refill helium gas of high purity (He 6.0, 99,9999%). Use the Service Kit. 	Bruker Service
35.040 hours (4 years)	Cryogenic Refrigerator	Repeat procedure after 17.520 hours (2 years).	Bruker Service
52560 hours (6 years)	Cryogenic Refrigerator	<ul style="list-style-type: none"> Replace rotary valve and cold head. Refer to chapter "Replacement of the Rotary Valve and the Cold Head" on page 7-118. Replace the compressor. Pump and flush the flex lines with helium gas of high purity (He 6.0, 99,9999%) 	Bruker Service

Table 7.2: Maintenance Timetable - part 2

7.4 Maintenance Work at the Cryogenic Refrigerator

Persons in charge: Bruker Service only

8 Disassembling

8.1 Safety

Persons in charge: Bruker Service only



A Appendix

Warning Signs

Figures

Tables

Glossary / Abbreviations

Technical Data

Index

A.1 Warning Signs

Gefahr/Danger

Art und Quelle der Gefahr	
Kind and Source of Danger	19

Warnung/Warning

Art und Quelle der Gefahr	
Kind and Source of Danger	20, 21, 22, 23, 24, 25, 26

Vorsicht/Caution

Art und Quelle der Gefahr	
Kind and Source of Danger	19, 27

Information

Hinweise, die dem Anwender die Arbeit erleichtern	
Information and links for trouble-free handling and operation	19

A.2 Figures

Figure 1.1:	General view Cryostat of the Magnet System.....	12
Figure 2.1:	Safety Devices of the Cryostat.....	28
Figure 3.1:	Packaging (without surrounding panels).....	31
Figure 3.2:	Transportation by forklift	33
Figure 3.3:	Attachment points for lifting equipment.....	34
Figure 5.1:	Start the Magnet Stand	40
Figure 5.2:	Stop the Magnet Stand	40
Figure 6.1:	Quench picture.....	53
Figure 6.2:	Temperature profile during Cryogenic Refrigerator failure.....	55
Figure A.1:	Dimension Cryostat - Front View	76
Figure A.2:	Dimension Cryostat - Top View	78
Figure A.3:	Helium Level Graph	80
Figure A.4:	Helium Level Sensor	81
Figure A.5:	Characteristic Curve PT 100.....	82
Figure A.6:	Characteristic Curve IBT	83
Figure A.7:	Wiring Diagramm Temperature Sensors	84
Figure A.8:	Current Lead 55 pins	86
Figure A.9:	Wiring Diagram Magnet	87
Figure A.10:	Wiring Diagram Magnet - Control	88
Figure A.11:	Wiring Diagram Magnet - Diagnostic and Temperature Sensors	89
Figure A.12:	Shorting Plug 55 pins	97
Figure A.13:	Fringe field plot	98

A.3 Tables

Table 7.1:	Maintenance Timetable - part 1	58
Table 7.2:	Maintenance Timetable - part 2	59
Table A.1:	Glossary	71
Table A.2:	Abbreviations	71
Table A.1:	Weight Cryostat.....	73
Table A.2:	Weight Noise Protection Cover	73
Table A.3:	Weight Cryogenic Refrigerator Parts	73
Table A.4:	Weight Noise Protection Cover	73
Table A.5:	Dimensions of Box for the Cryostat	74
Table A.6:	Dimensions of Box for the Cryogenic Refrigerator.....	74
Table A.7:	Dimensions of Box for the Noise Protection Cover	74
Table A.8:	Environmental conditions	75
Table A.9:	Dimension Cryostat - Front View	77
Table A.10:	Dimensions Cryostat - Top View	78
Table A.11:	Installation Consumption.....	79
Table A.12:	Cryogenic Agents.....	79
Table A.13:	Helium Level Sensor	81
Table A.14:	Characteristic Values PT 100.....	82
Table A.15:	Characteristic Values IBT	83
Table A.16:	Specification of the Magnet.....	85
Table A.17:	Resistance at Room Temperature	91
Table A.18:	Heater currents	92
Table A.19:	Energizing assignment and currents.....	92
Table A.20:	Energizing currents	93
Table A.21:	Cycling assignment (recommended only for 500Mhz and more).....	94
Table A.22:	Deenergizing Assignment and Currents	95
Table A.23:	Fringe Field Values	98

A.4 Glossary / Abbreviations

Used term	Description
Box	Any kind of package used to protect sensitive parts during transportation.
Cryostat	The collective of all parts providing a temperature of 4 K inside for the superconducting magnet. The cryostat also provides the safety devices and the access ports for the cryogenic agents and electricity. The superconducting magnet inside the cryostat is not energized.
Dewar	Any kind of package used for transporting cryogenic agents like gaseous or liquid helium or nitrogen.
Magnet System	The collective of all parts necessary for the intended use. The superconducting magnet inside the cryostat is energized.

Table A.1: Glossary

Abbreviations	Description
ACD	Automatic Cooling Device
BSMS	Bruker Smart Magnetcontrol System
BSVT	Bruker Smart Variable Temperature System
MICS	Magnet Information and Control System; Bruker BioSpin registered trademark
RS	Radiation Shield
RT	Room Temperature; used for room temperature level and as prefix of parts witch are at room temperature
VTA	Variable Temperature Adapter
USR	Ultra Shielded Refrigerated; Bruker BioSpin registered trademark

Table A.2: Abbreviations

A.5 Technical Data MS 400'54 USR

A.5.1 Dimensions

Weights

Cryostat	Value	Unit
Weight, empty, without magnet stand	386	kg
Weight, completely filled, without magnet stand	405	kg
Weight, empty, ready for transportation including box	535	kg

Table A.1: Weight Cryostat

Magnet Stand	Value	Unit
Weight, ready for use	128	kg
Weight, ready for transportation including box	170	kg

Table A.2: Weight Noise Protection Cover

Cryogenic Refrigerator Parts	Value	Unit
Weight, Cryogenic Refrigerator Parts	80	kg
Weight, Rotary Valve Column	65	kg
Weight, ready for transportation including box	220	kg

Table A.3: Weight Cryogenic Refrigerator Parts

Noise Protection Cover	Value	Unit
Weight, ready for use	25	kg
Weight, empty, ready for transportation including box	46	kg

Table A.4: Weight Noise Protection Cover

Dimensions of Transportation Boxes

Cryostat	Value	Unit
Height	1800	mm
Width	1260	mm
Depth	1050	mm

Table A.5: Dimensions of Box for the Cryostat

Cryogenic Refrigerator Parts	Value	Unit
Height	950	mm
Width	790	mm
Depth	1150	mm

Table A.6: Dimensions of Box for the Cryogenic Refrigerator

Noise Protection Cover	Value	Unit
Height	940	mm
Width	800	mm
Depth	1200	mm

Table A.7: Dimensions of Box for the Noise Protection Cover

Environmental conditions

	Value	Unit
Minimum surrounding temperature	7	°C
Maximum surrounding temperature	38	°C
Maximum relative humidity up to 31°C	80	%
Maximum relative humidity between 31°C and 40°C linear decreasing	80-50	%

Table A.8: Environmental conditions

Identification Plate

The identification plate is on the right back side of the cryostat fixed at the bottom plate.

Contents of the identification plate:

- Address of the Manufacturer
- Magnet System Identifier
- Type
- Identification Number
- Magnet Identifier
- Serial Number
- Year of Construction
- Cryostat Identifier
- Specification Helium Vessel
- Specification Vacuum Chamber
- Weight (empty and completely filled) including magnet stand

Dimensions Cryostat

Front view

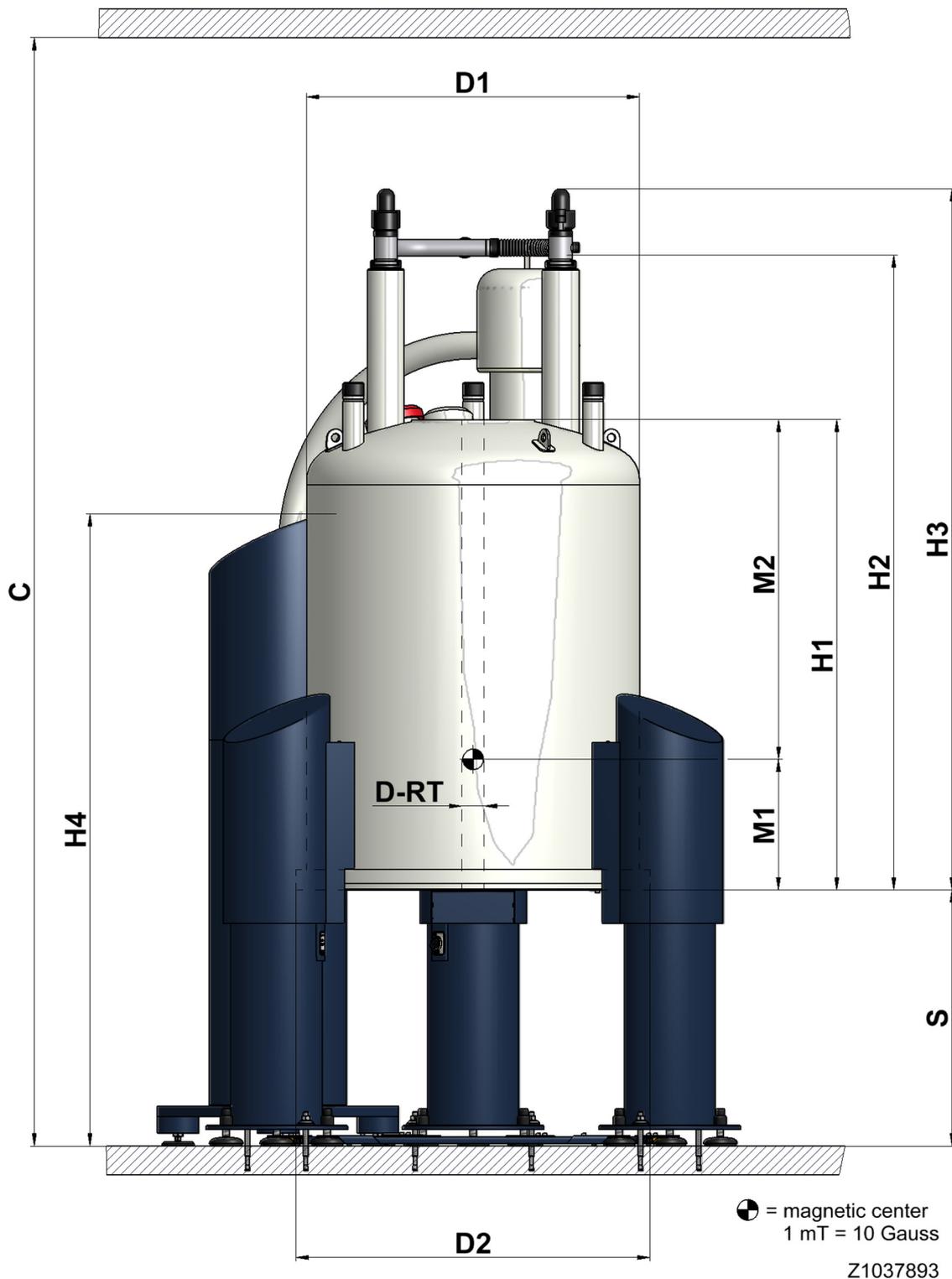


Figure A.1: Dimension Cryostat - Front View

Dimensions Cryostat	Value	Unit
C Operational Ceiling Height	3110	mm
D-RT Diameter RT Bore Tube	54	mm
D1 Diameter RT vessel	800	mm
D2 Diameter Bottom Plate	850	mm
H1 Height Cryostat (bottom plate to top flange)	1140	mm
H2 Height Cryostat (minimum height for transportation)	1538	mm
H3 Height Cryostat (bottom plate to flow system)	1700	mm
H4 Height Rotary Valve Column	1530	mm
S Height Magnet Stand (floor to bottom plate)	620	mm
M1 Distance magnetic Center to bottom flange (calculated)	317	mm
M2 Distance magnetic center to top flange (calculated)	823	mm

Table A.9: Dimension Cryostat - Front View

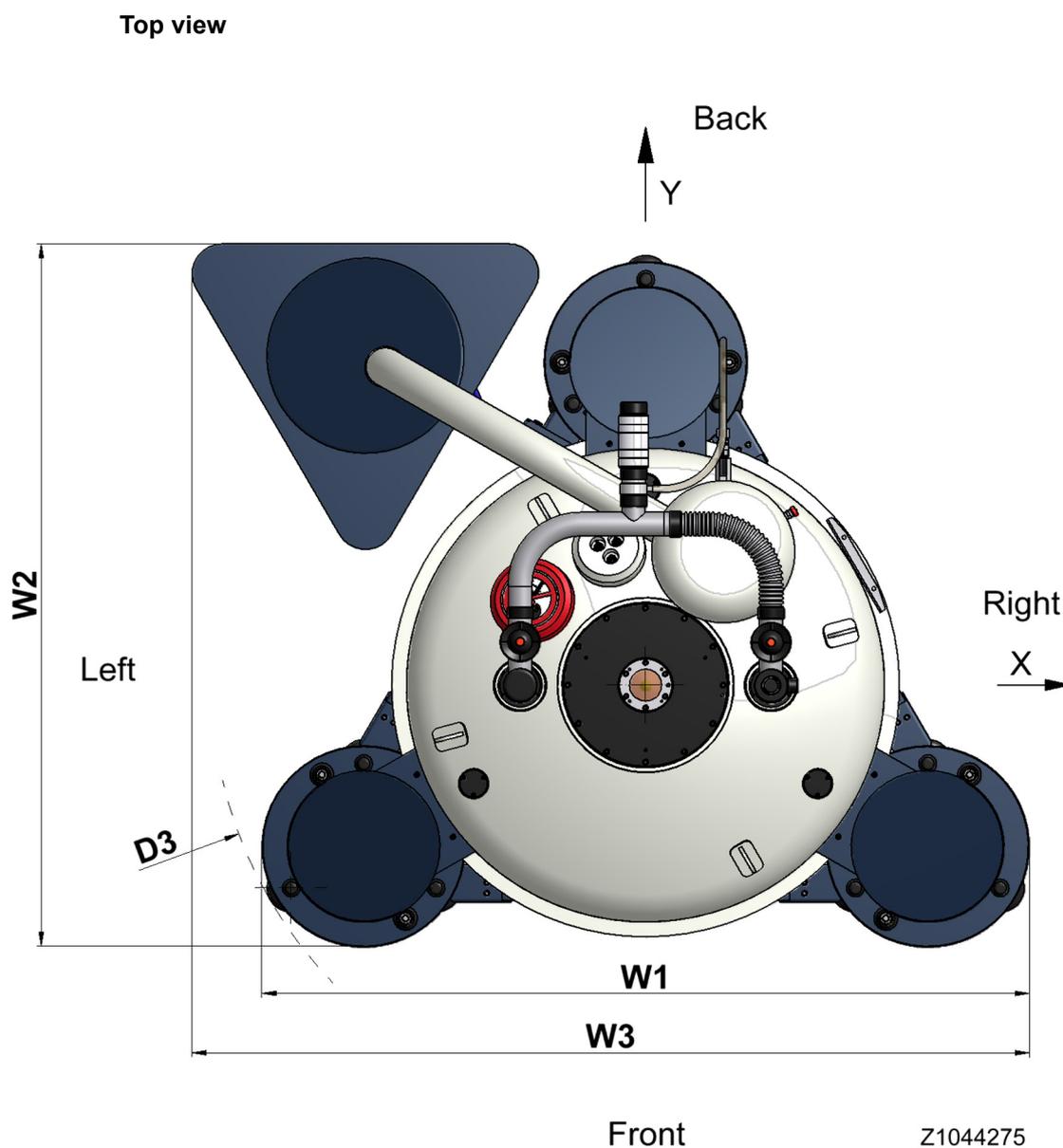


Figure A.2: Dimension Cryostat - Top View

Dimensions Cryostat	Value	Unit
W1	1284	mm
W2	1186	mm
W3	1401	mm
D3 ¹	1451	mm

Table A.10: Dimensions Cryostat - Top View

1. Keep at least an additional free space of 1.5 m around the magnet system for service.

A.5.2 Technical Data Cryostat

Supplies during installation

The consumption of liquid cryogenic agents during installation consists of consumption for cooling down the cryostat and the volume to fill the vessels.

Cryogenic Agent	Value	Unit
Nitrogen	500	l
Helium	400	l

Table A.11: Installation Consumption

Fill Volume, Evaporation Rate and Hold Time

Cryogenic Agents	Value	Unit
Helium vessel total volume	148	l
Helium refill volume	116	l
Helium evaporation rate	13	ml/h
Helium hold time ¹ with Cryogenic Refrigerator operating	372	days
Helium hold time without Cryogenic Refrigerator operating (at minimum allowed helium level)	3	days
Helium refill volume after Quench (cool down and refill)	150	l
(Maximum allowed pressure difference) * (Volume of the helium vessel) ²	0.6*148 = 89	bar*l

Table A.12: Cryogenic Agents

1. Maximum time intervall between two fillings

2. Maximum allowed operation pressure 200 mbar (4.9 psi), maximum allowed pressure between the vessel and the RT vacuum chamber 1200 mbar (19 psi)

Helium Level Graph

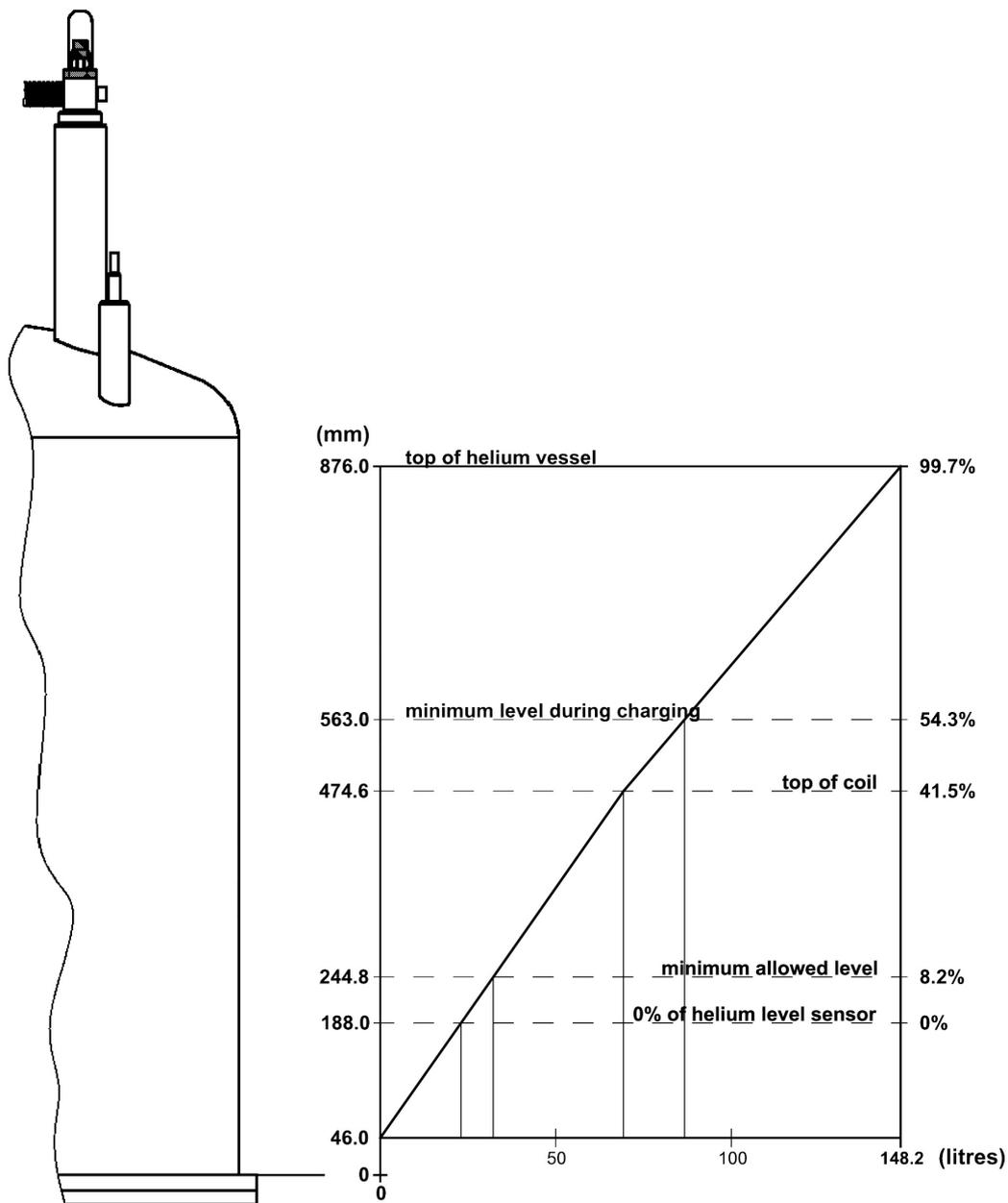



Figure A.3: Helium Level Graph

Helium Level Sensor

The helium level sensor is inserted in the helium fill-in turret.

Helium Level Sensor	Material No.	Value	Unit
Level Sensor Type	Z58114	1300/690	
Overall length		1300	mm
Active length		690	mm
Calibration 0 %, Calibration resistor	Z28630 blue	140	Ω
Calibration 100 %, Calibration resistor	Z28628 black	15	Ω

Table A.13: Helium Level Sensor

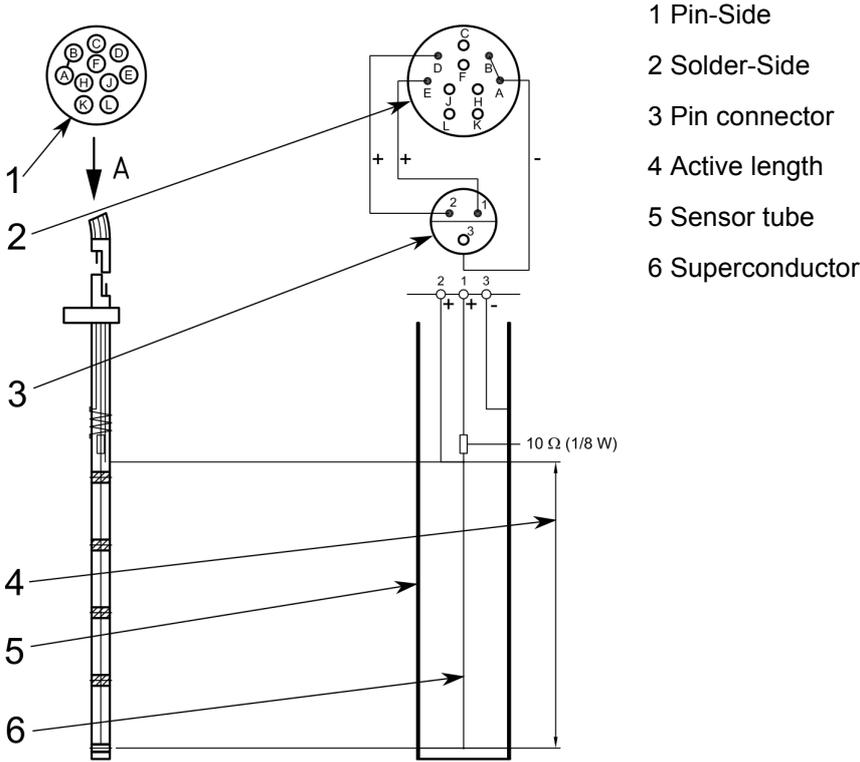

Z1029564

Figure A.4: Helium Level Sensor

Temperature Sensors

The temperature sensors (PT 100 and IBT) will be used to monitor the temperature of the magnets during cooling and warming up of the magnet system.

PT 100

i Measure the resistance with a maximum current of 1 mA.

	Temperature	Unit	Resistance	Unit
Room Temperature	293	K	107.8	Ω
	273	K	100.0	Ω
	250	K	91.0	Ω
	200	K	71.1	Ω
	150	K	50.9	Ω
	100	K	30.0	Ω
Liquid Nitrogen	77	K	20.1	Ω

Table A.14: Characteristic Values PT 100

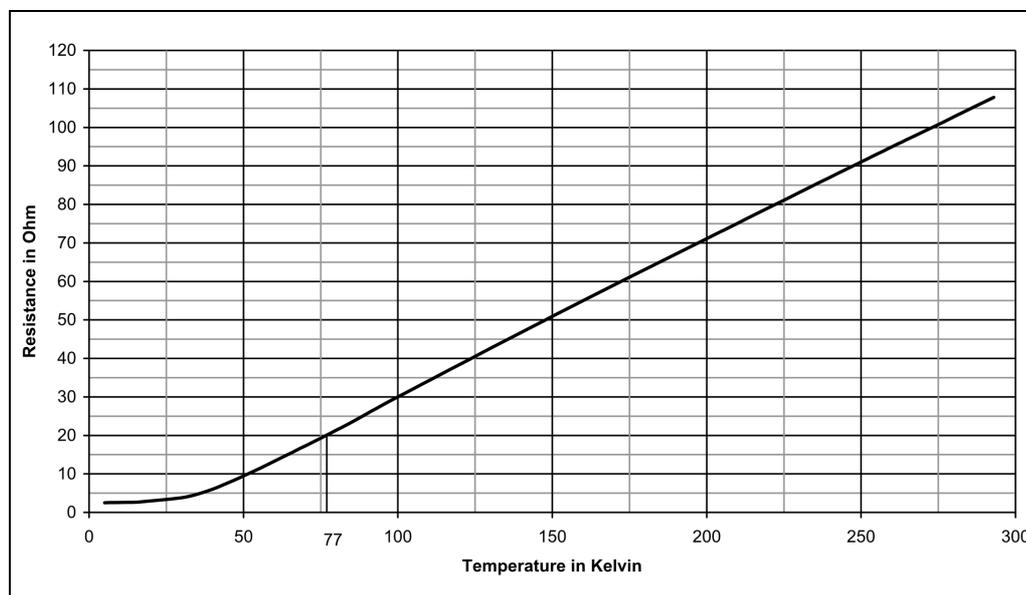


Figure A.5: Characteristic Curve PT 100

IBT Carbon Resistor

i Measure the Resistance with a maximum current of 0.1 mA.

	Temperature	Unit	Resistance	Unit
Room Temperature	300	K	220	Ω
Liquid Nitrogen	77	K	265	Ω
	40	K	300	Ω
	20	K	350	Ω
	10	K	420	Ω
	8	K	450	Ω
	6	K	500	Ω
	5	K	540	Ω
Liquid Helium	4.2	K	575	Ω

Table A.15: Characteristic Values IBT

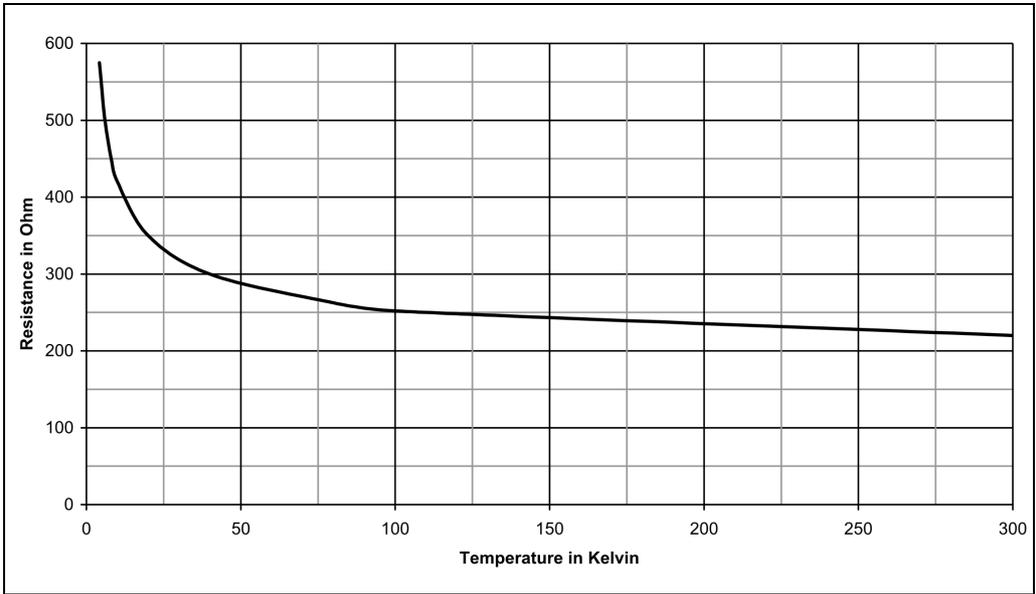


Figure A.6: Characteristic Curve IBT

Wiring Diagramm Temperature Sensors

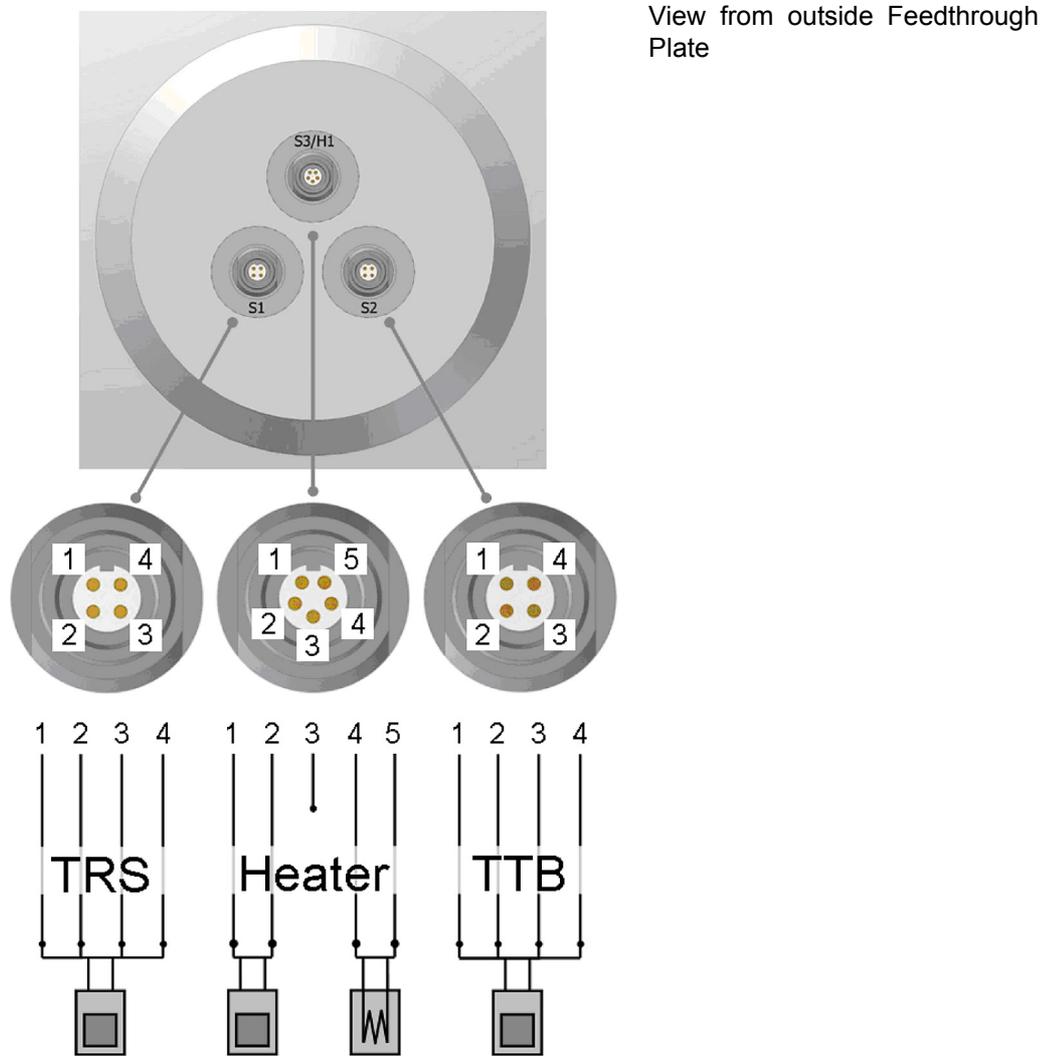


Figure A.7: Wiring Diagramm Temperature Sensors

A.5.3 Technical Data Magnet

Technical Data Magnet	Value	Unit
Proton frequency	400	MHz
Central field	9.4	T
Coil inductance	39.6	H
Magnetic energy	210	kJ
Magnetic centre from bottom flange (design value)	317	mm
Maximum drift rate	0.01 4	ppm/h Hz/h
Radial stray field at 0.5mT	0.5	m
Axial stray field at 0.5mT	1	m

Table A.16: Specification of the Magnet

Operating modes of the Magnet System:


Driven Mode

In the driven mode the current lead is mounted and the electricity is flowing through the power supply. The coils of the magnet can be charged and discharged.

Persistent Mode

In the persistent mode the electricity is flowing exclusively within the magnet. The circuit has no connection to the outside. The magnetic field cannot be switched off.


Current Lead

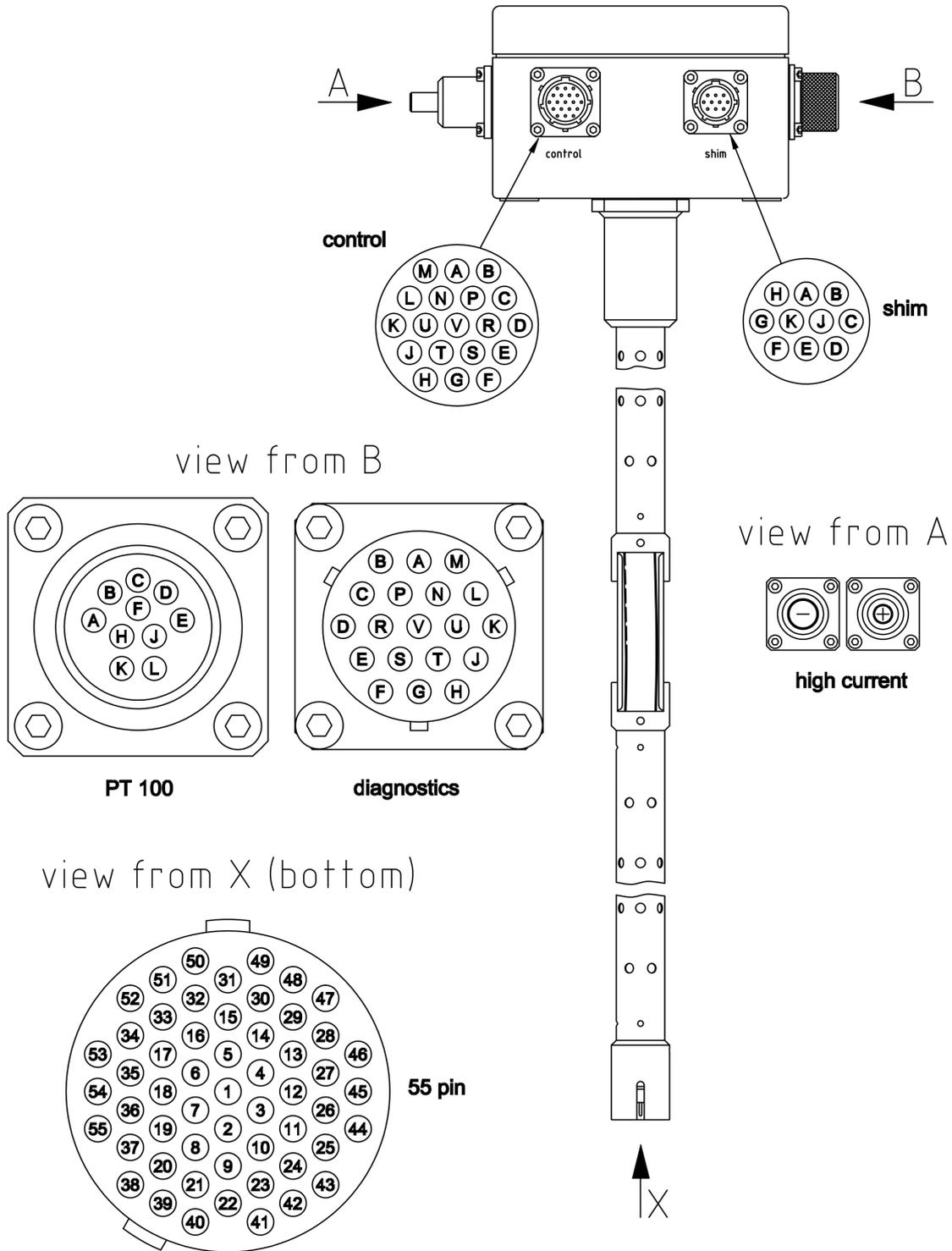

Z1029367

Figure A.8: Current Lead 55 pins


Wiring Diagram Magnet

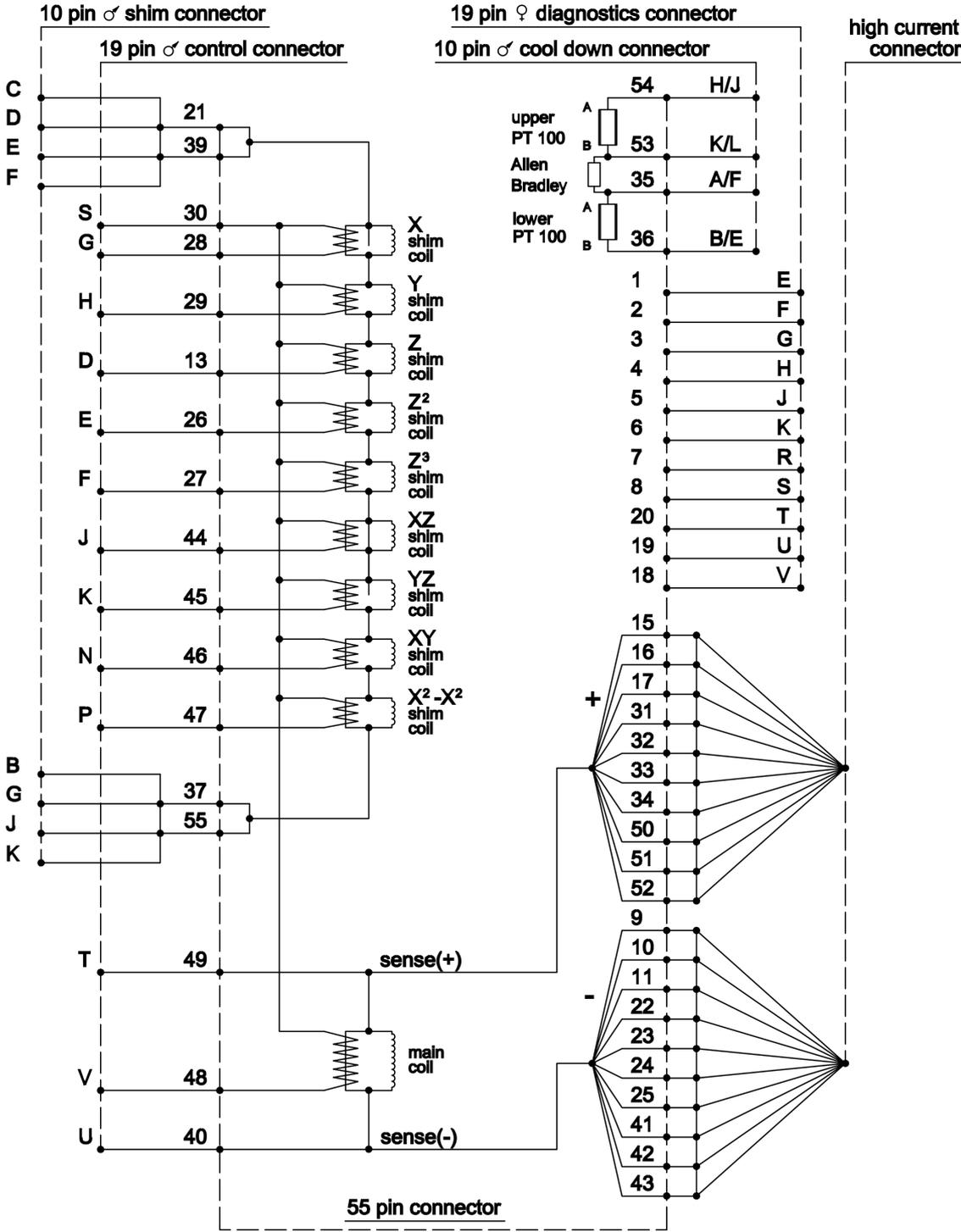

Z1030082

Figure A.9: Wiring Diagram Magnet


Wiring Diagram Magnet Control

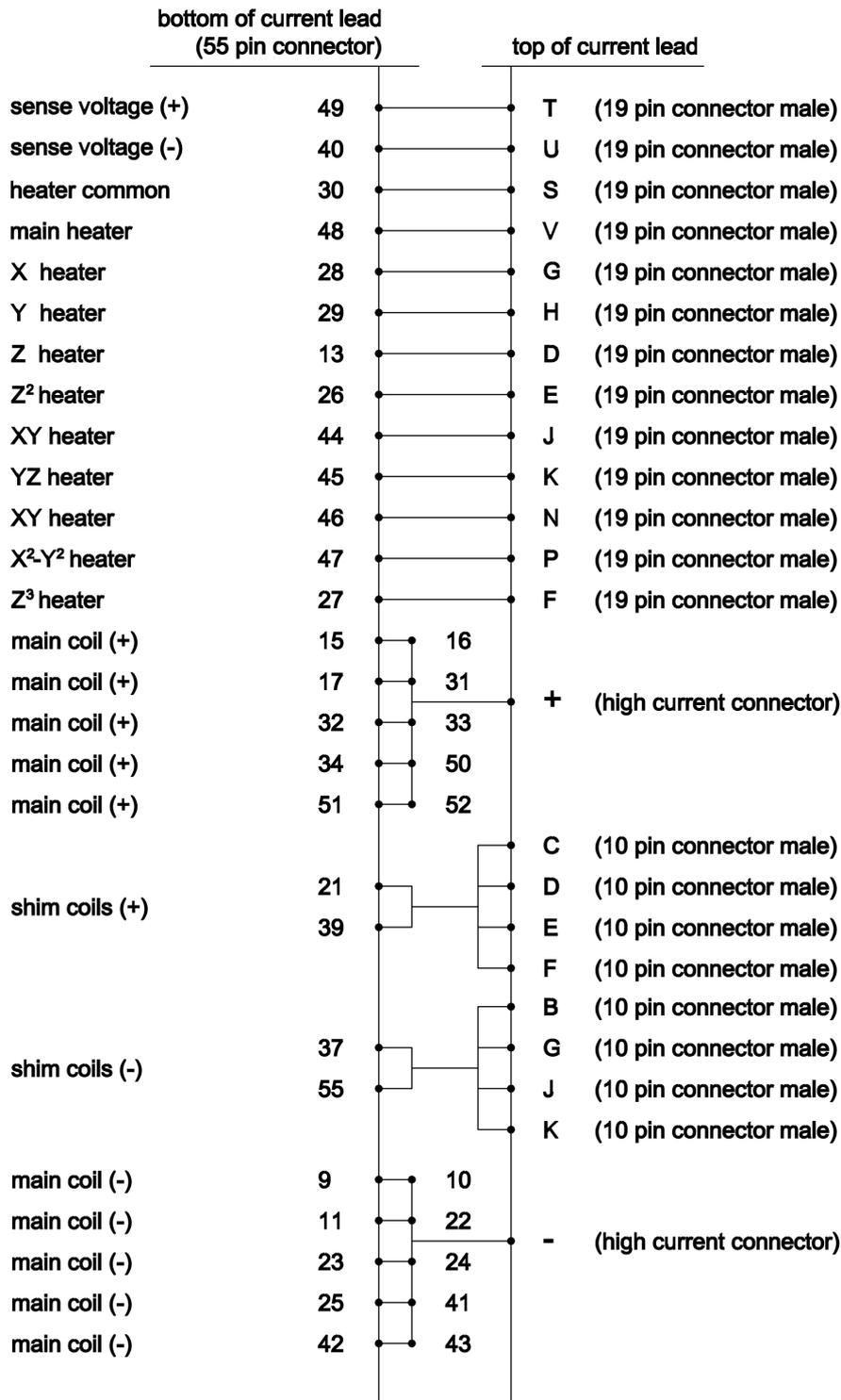

Z1029391

Figure A.10: Wiring Diagram Magnet - Control

Wiring Diagram Magnet Diagnostic and Temperature Sensors

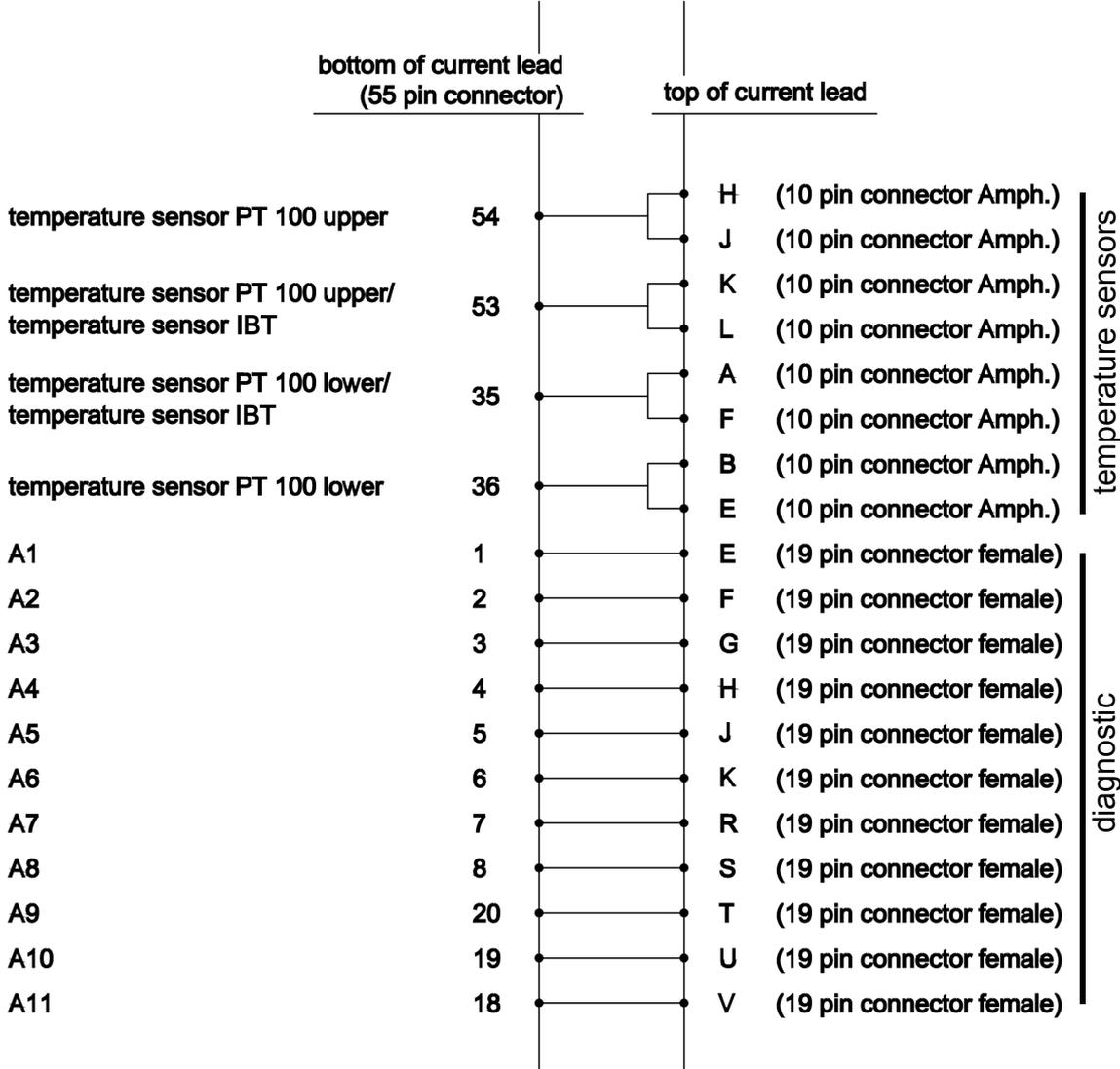

Z1029392

Figure A.11: Wiring Diagram Magnet - Diagnostic and Temperature Sensors

This side is intentionally left blank.

Resistance at Room Temperature

	Pin	Connector	Description	Value	Unit
From: To:	PIN V PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	Main Heater		Ω
From: To:	PIN D PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	Z Heater		Ω
From: To:	PIN G PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	X Heater		Ω
From: To:	PIN H PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	Y Heater		Ω
From: To:	PIN J PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	XZ Heater		Ω
From: To:	PIN K PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	YZ Heater		Ω
From: To:	PIN N PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	XY Heater		Ω
From: To:	PIN P PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	X ² -Y ² Heater		Ω
From: To:	PIN E PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	Z ² Heater		Ω
From: To:	PIN F PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	Z ³ Heater		Ω
From: To:	PIN C,D,E,F PIN B,G,J,K	10 PIN Con SHIM 10 PIN Con SHIM	Shim Coils +/-		Ω
From: To:	+ PIN T	High current Con 19 PIN Con CONTROL	High Current to Sense +		Ω
From: To:	+ -	High current Con High current Con	Main Coil		Ω
From: To:	- PIN U	High current Con 19 PIN Con CONTROL	High Current to Sense -		Ω
From: To:	PIN T PIN U	19 PIN Con CONTROL 19 PIN Con CONTROL	Sense + Sense -		Ω
From: To:	PIN C PIN S	10 PIN Con SHIM 19 PIN Con CONTROL	Shim Coil to Heater (common)		Ω
From: To:	PIN C PIN T	10 PIN Con SHIM 19 PIN Con CONTROL	Shim Coil to Main Coil		Ω
From: To:	PIN T PIN S	19 PIN Con CONTROL 19 PIN Con CONTROL	Sense to Heater (common)		Ω
From: To:		All Connectors Ground	Insulation Magnet to Cryostat		Ω
From: To:	PIN K PIN H	10 Pin cool down Con 10 Pin cool down Con	Upper temperature sensor PT 100		Ω
From: To:	PIN A PIN K	10 Pin cool down Con 10 Pin cool down Con	IBT Carbon temperature sensor		Ω
From: To:	PIN A PIN B	10 Pin cool down Con 10 Pin cool down Con	Lower temperature sensor PT100		Ω

Table A.17: Resistance at Room Temperature

Heater Currents



Heater Currents	Value	Unit
Main heater current		A
Shim heater current		A

Table A.18: Heater currents

Energizing Assignment and Currents

Energizing Currents [A]			Sense voltage [mV]	Remarks Bruker Test Site
0	to	120	4000	
120	to	170	2000	
170	to	190	1000	
190	to	205	500	
205	to		200	
Overshoot (0.25 % of final current)				
	to		50	
10 Minutes break at overshoot current				
	to		-50	
Total energizing time [min]				

Table A.19: Energizing assignment and currents

Energizing Currents



Energizing Currents	Value Bruker Test Site	Value Customer Site #1	Value Customer Site #2	Unit
Magnet main current				A
Z-Shim current				A
Z ² -Shim current				A
Z ³ -Shim current				A
X-Shim current				A
Y-Shim current				A
XZ-Shim current				A
YZ-Shim current				A
XY-Shim current				A
X ² -Y ² -Shim current				A
Magnetic center from bottom flange (measured)				mm
Magnetic center from top flange (measured)				mm
Date and Signature				

Table A.20: Energizing currents

Cycling Assignment and Shim Currents

(recommended only for 500MHz and more)



Cycling Assignment and Shim Currents	Value	Unit
Time between energizing and cycling (at least 12 h)		h
Time between cycling and shimming (at least 12 h)		h
Current rate		A/h
Z-Shim current		A
Z ² -Shim current		A
Z ³ -Shim current		A
X-Shim current		A
Y-Shim current		A
XZ-Shim current		A
YZ-Shim current		A
XY-Shim current		A
X ² -Y ² -Shim current		A
Frequency change due to cycling		kHz
Date and Signature		

Table A.21: Cycling assignment (recommended only for 500Mhz and more)

Deenergizing Assignment and Currents




Deenergizing Currents [A]			Sense voltage [mV]	Remarks Bruker Test Site
	to	205	400	
205	to	190	1000	
190	to	0	2000	through external diodes
Total deenergizing time [min]				

Table A.22: Deenergizing Assignment and Currents

This side is intentionally left blank.

Shorting Plug

The shorting plug is plugged after removal of the current lead. After inserting the shorting plug the current flows through the shorting plug and no longer through the current lead and the power supply.

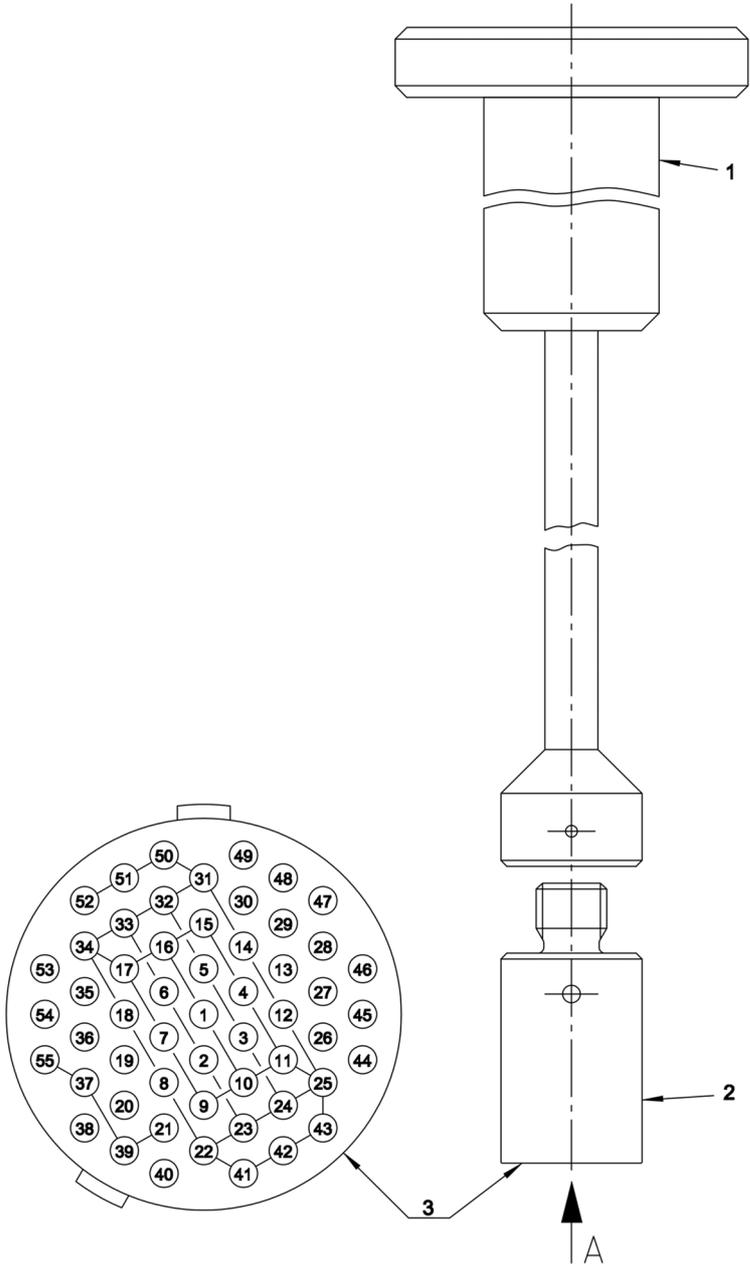

Z1029521

Figure A.12: Shorting Plug 55 pins
1 Shorting Plug Tool for fitting and removing the shorting plug
2 Shorting Plug
3 Shorting Plug - view from pin side

Fringe Field Plot

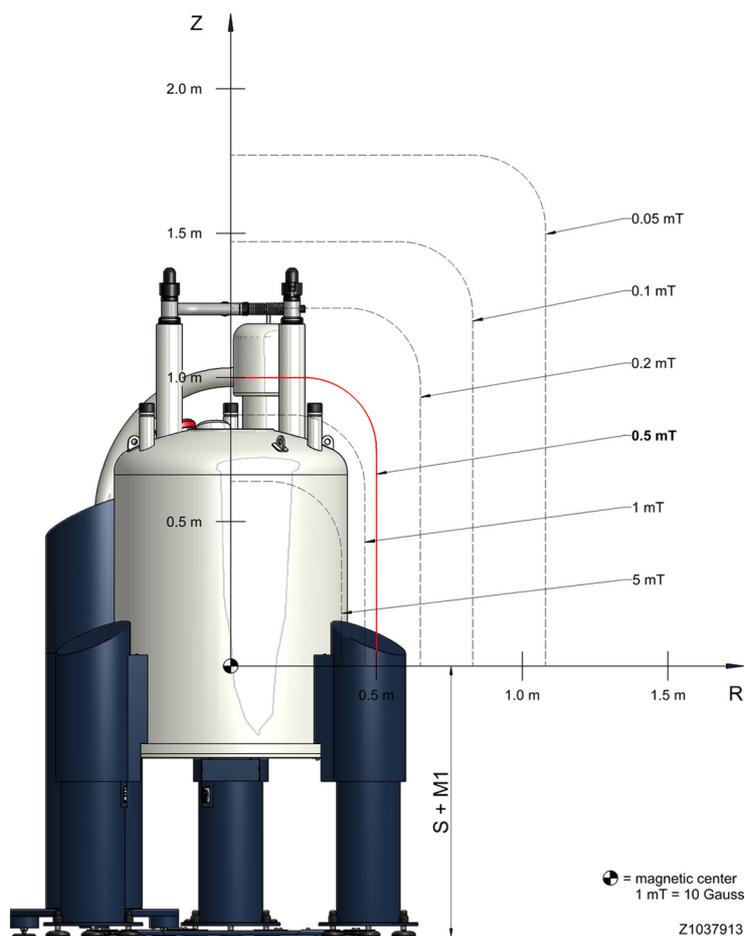


Figure A.13: Fringe field plot

Fringe Field	Unit	R max	Unit	Z max	Unit
0.05	mT	1.08	m	1.77	m
0.1	mT	0.83	m	1.47	m
0.2	mT	0.65	m	1.24	m
0.5	mT	0.50	m	1.00	m
1.0	mT	0.46	m	0.87	m
3.0	mT	0.41	m	0.70	m
5.0	mT	0.38	m	0.64	m

Table A.23: Fringe Field Values

A.6 Index

A

Approved Persons of the Customer	15
Attachment Points	34

B

Bruker Service	15
----------------------	----

C

Cleaning Staff	16
Cold Head	12
Complaint Process	32
Compressor	12
Connection Line	12
Cryogenic Agents	20, 21
Cryogenic Refrigerator	13, 55
Customer Responsibilities	16

D

Disturbances	43
Drop Off Plate	28

E

Electricity	21
Emergency	29
Emergency Exit	27

F

Fire Department	29
First Aid for Cold Burns	29
Flex Lines	12

I

Identification Plate	75
Intended Use	18, 26

L

Liability	11
Low Temperatures	23

M

Magnet Stand	12, 40
Magnetic Field	20

P

Portable Oxygen Monitor	26
Protective Clothes	26
Protective Gloves	26
Protective Goggles	26

Q

Quench	22, 53
Quench Valve	28

R

Rescue Equipment	29
Rotary Valve Column	12

S

Safety Devices	19
Safety Shoes	26
Shock Watch	32
Slippage	23
Spontaneous Ignition	23

T

Tilt Watch	32
Tilting	24

W

Warranty	11
----------------	----

Revision History List

Index:	Date:	Alteration Type:
00	February 2008	First release, design of the magent system not yet fixed.
01	April 2010	New release after re-design of the cryostat and release of the Rotary Valve Column, using a new specific form of the Corporate Guide Line and the definition from the Bruker documentation group.
02	November 2010	New release with a new order of content.



Bruker BioSpin

your solution partner

Bruker BioSpin provides a world class, market-leader range of analytical solutions for your scientific and materials science needs. Our solution-based approach enables us to work closely with you to identify your specific demand and determining the best solution package from our comprehensive range. Furthermore, we offer you the exclusive opportunity to collaborate with us on new developments.

Our ongoing efforts and substantial investment in research and development demonstrate our long-term commitment to technological innovation on behalf of our customers. With 50 years of experience in meeting the demands of the scientific community across a range of specialist disciplines, Bruker BioSpin has built a sound rapport with professionals from the scientific community through understanding their specific demands and providing them with effective solutions and a service of unmatched quality.

● **Bruker BioSpin Group**

info@bruker-biospin.com
www.bruker-biospin.com